R453Plus1Toolbox

March 24, 2012

AVASet-class

Class to contain Amplicon Variant Analyzer Output

Description

Container to store data imported from a project of Roche's Amplicon Variant Analyzer Software. It stores all information into an extended version of the Biobase ExpressionSet.

Objects from the Class

Objects can be created by calls of the form AVASet (filename). While filename is the path of the project data (i.e. a directory that contains the files and subdirectories "Amplicons/ProjectDef/ampliconsProject.txt", "Amplicons/Results/Variants/currentVariantDefs.txt", "Amplicons/Results/Variants", "Amplicons/Results/Align").

Slots

- assayData: Object of class AssayData. Contains the number of reads and the total read depth for every variant and each sample in forward and reverse direction. Its column number equals nrow (phenoData).
- featureData: Object of class AnnotatedDataFrame. Contains information about the type, position and reference of each variant.
- phenoData: Object of class AnnotatedDataFrame. Contains the sample-IDs and name, annotation and group of the read data for all samples. If available, the lane, pico titer plate(s) (PTP) or MID(s) of each sample are shown as well.
- assayDataAmp: Object of class AssayData. Contains the number of reads for every amplicon and each sample in forward/reverse direction. Its column number equals nrow (featureDataAmp).
- featureDataAmp: Object of class AnnotatedDataFrame. Contains the primer sequences, reference sequences and the coordinates of the target regions for every amplicon.
- referenceSequences: Object of class AlignedRead. If additional alignment information were computed via alignShortReads, this slot knows about the chromosome, position and the strand of each reference sequence.
- variantFilterPerc: Object of class numeric. Contains a threshold to display only those variants, whose coverage (in percent) in forward and reverse direction in at least one sample is higher than this filter value. See setVariantFilter for details about setting this value.
- variantFilter: Object of class character. Contains a vector of variant names whose coverage (in percent) in forward and reverse direction in at least one sample is higher than the filter value in variantFilterPerc.

2 AVASet-class

dirs: Object of class character. Based on a directory given at instantiation of the object, it contains a vector of several directories containing all relevant AVA-project files.

- experimentData: Object of class MIAME. Contains details of the experiment.
- annotation: Object of class character. Label associated with the annotation package used in the experiment.
- protocolData: Object of class annotatedDataFrame. Contains additional information about the samples.
- .__classVersion__: Object of class Versions. Remembers the R and R453Toolbox version numbers used to created the AVASet instance.

Extends

Class eSet, directly. Class VersionedBiobase, by class "eSet", distance 2. Class Versioned, by class "eSet", distance 3.

Methods

- **object[i,j**:] Allows subsetting an AVASet object by features (i) and samples (j).
- **assayDataAmp(object), assayDataAmp(object)<-value:** Similar to assayData of the Biobase ExpressionSet, this function returns/replaces the amplicon assay data.
- **fDataAmp(object):** Similar to fData of the Biobase ExpressionSet, this function returns the amplicon feature data as a data frame.
- **featureDataAmp(object), featureDataAmp(object)<-value:** Similar to featureData of the Biobase ExpressionSet, this function returns/replaces the amplicon feature data and feature meta.
- **referenceSequences(object), referenceSequences(object)<-value:** Returns/replaces the reference sequence slot.
- **alignShortReads(object, bsGenome):** Retrieve the chromosomal positions of the amplicon sequences.
- **setVariantFilter(object):** Sets the filter to display only those variants, whose coverage (in percent) in forward and reverse direction in at least one sample is higher than the given value.
- **getVariantPercentages(object)** Computes the coverage for every variant over all reads (forward and/or reverse) and for each sample.
- annotateVariants(object): Annotates given genomic variants. See annotateVariants for details.
- htmlReport(object): Exports all (filtered) variant data into a html report. See htmlReport for details

Author(s)

Christoph Bartenhagen

See Also

MapperSet-class, annotateVariants, alignShortReads, htmlReport, setVariantFilter, getVariantPercentages

AVASet 3

Examples

```
# sum up class structure
showClass("AVASet")

# load an AVA dataset containing 6 samples, 4 amplicons and 259 variants
data(avaSetExample)
avaSetExample

# show contents of assay, feature and pheno data
head(assayData(avaSetExample)$variantForwCount)
head(assayData(avaSetExample)$totalForwCount)
head(assayData(avaSetExample)$variantRevCount)
head(assayData(avaSetExample)$totalRevCount)
head(fData(avaSetExample))
pData(avaSetExample)
assayDataAmp(avaSetExample)
fDataAmp(avaSetExample)
referenceSequences(avaSetExample)
```

AVASet

Creating an AVASet

Description

This function imports a project of Roche's Amplicon Variant Analyzer Software. It stores all information into an extended version of the Biobase eSet.

Usage

AVASet (dirname)

Arguments

dirname

The path of the project data (i.e. a directory that contains the files and subdirectories "Amplicons/ProjectDef/ampliconsProject.txt", "Amplicons/Results/Variants/currentVariantDef "Amplicons/Results/Variants", "Amplicons/Results/Align").

Details

An AVASet object consists of three slots to store data about 1. variants

variantForwCount/variantRevCount: Data frames that contain the number of reads with the respective variant in forward/reverse direction.

totalForwCount/totalRevCount: Data frames that contain the total coverage for every variant location in forward/reverse direction.

referenceSeq: Gives the identifier of the reference sequence.

variantBase/referenceBases: The bases changed in each variant.

4 AVASet

start/end: The position of the variant on the reference sequence.

canonicalPattern/name: Short identifiers of a variant including the position and the bases changed.

2. amplicons

forwCount/revCount: Data frames that contain the number of reads for every amplicon and each sample in forward/reverse direction.

primer1,primer2: The primer sequences for every amplicon.

referenceSeqID: The identifier of the reference sequence.

targetStart/targetEnd: The coordinates of the target region.

3. reference sequences

If additional information has been loaded from Ensembl via alignShortReads, this slot knows about the chro

The structure of the variant and amplicon data is derived from the Biobase eSet and thus separated into assayData, phenoData and featureData. All information about the reference sequences is stored into an object of class AlignedRead.

The phenoData of the variants lists the sample-IDs and name, annotation and group of the read data for all samples. If available, the pico titer plate(s) (PTP) or MID(s) of each sample are shown as well.

Value

An instance of the AVASet class.

Author(s)

Christoph Bartenhagen

See Also

MapperSet-class, alignShortReads

Examples

```
# load an AVA dataset containing 6 samples, 4 amplicons and 259 variants
data(avaSetExample)
avaSetExample
```

AnnotatedVariants-class 5

AnnotatedVariants-class

Class "AnnotatedVariants"

Description

A class for storing annotation about variants. An object of this class is returned by the method annotateVariants. The class has not been designed to be created by users directly.

Details

The list encapsulated by this class has one element for each variant. Each element is a nested list with the elements genes, transcripts, exons and snps. All these elements are data frames listing genes, transcripts, exons or snps respectively that were affected by the variant. Use the example below to explore the data frames' contents.

Objects from the Class

Objects can be created by calls of the form new("AnnotatedVariants"). The method annotateVariants returns AnnotatedVariants-objects.

Slots

annotated Variants: Object of class "list" with one entry for each variant.

Methods

```
annotatedVariants signature(object = "AnnotatedVariants"): Get the list with
    variants.

annotatedVariants<- signature(object = "AnnotatedVariants", value = "list"):
    Set a new list with variants.

names signature(x = "AnnotatedVariants"): Get the names of the with variants.

names<- signature(x = "AnnotatedVariants", value = "character"): Set the
    names of the variants.</pre>
```

Author(s)

Hans-Ulrich Klein

See Also

annotateVariants, htmlReport

Examples

```
variants = data.frame(
    start=c(106157528, 106154991,106156184),
    end=c(106157528, 106154994,106156185),
    chromosome=c("4", "4", "4"),
    strand=c("+", "+", "+"),
    seqRef=c("A", "ATAG", "---"),
    seqMut=c("G", "----", "ATA"),
```

6 Breakpoints-class

```
seqSur=c("TACAGAA", "TTTATAGATA", "AGC---TCC"),
    stringsAsFactors=FALSE)
rownames(variants) = c("snp", "del", "ins")
## Not run: av = annotateVariants(variants)
## Not run: annotatedVariants(av)[["snp"]]
```

```
Breakpoints-class Class "Breakpoints"
```

Description

Container to store chimeric reads that were clustered to putative breakpoints indicating structural variants. Related information like breakpoint position or alignment information about the chimeric reads is stored as well.

Objects from the Class

Objects can be created by calls of the form new("Breakpoints", ...). Usually, objects will be created by calling the detectBreakpoints method. It is not intended that users create objects of this class manually.

All slots of this class can be found twice. One slot name ends with "C1" and the other "C2". The slots labeled with "C2" are empty until mergeBreakpoints has been called and contain information about putatively associated breakpoints detected by mergeBreakpoints.

Slots

- seqsC1: Object of class "list" with one data frame for each breakpoint. The data frame stores all chimeric reads covering the first breakpoint together with the alignment information.
- seqsC2: Object of class "list" with one data frame for each breakpoint. The data frame stores all chimeric reads covering the second breakpoint together with the alignment information.
- commonBpsC1: Object of class "list" with one data frame for each breakpoint. The data frame stores the consensus breakpoint sequence as well as the breakpoint coordinates of the first breakpoint.
- commonBpsC2: Object of class "list" with one data frame for each breakpoint. The data frame stores the consensus breakpoint sequence as well as the breakpoint coordinates of the second breakpoint.
- commonAlignC1: Object of class "list" with one object of class PairwiseAlignedFixedSubject-class for each breakpoint storing the alignments of the chimeric reads against the consensus breakpoint sequence for the first breakpoint.
- commonAlignC2: Object of class "list" with one object of class PairwiseAlignedFixedSubject-class for each breakpoint storing the alignments of the chimeric reads against the consensus breakpoint sequence for the second breakpoint.
- alignedReadsC1: Object of class "list" with one object of class AlignedRead-class storing all chimeric reads covering the first breakpoint and their alignments.
- alignedReadsC2: Object of class "list" with one object of class AlignedRead-class storing all chimeric reads covering the second breakpoint and their alignments.

Breakpoints-class 7

Methods

```
alignedReadsC1<- signature(object = "Breakpoints", value = "list"): Setter-
method for the alignedReadsC1 slot.</pre>
```

- alignedReadsC1 signature(object = "Breakpoints"): Getter-method for the alignedReadsC1 slot.
- alignedReadsC2<- signature(object = "Breakpoints", value = "list"): Settermethod for the alignedReadsC2 slot.</pre>
- alignedReadsC2 signature(object = "Breakpoints"): Getter-method for the alignedReadsC2 slot.
- commonAlignC1<- signature(object = "Breakpoints", value = "list"): Settermethod for the commonAlignC1 slot.
- commonAlignC1 signature(object = "Breakpoints"): Getter-method for the commonAlignC1 slot.
- commonAlignC2<- signature(object = "Breakpoints", value = "list"): Settermethod for the commonAlignC2 slot.
- commonAlignC2 signature(object = "Breakpoints"): Getter-method for the commonAlignC2 slot.
- commonBpsC1<- signature(object = "Breakpoints", value = "list"): Settermethod for the commonBpsC1 slot.</pre>
- commonBpsC2<- signature(object = "Breakpoints", value = "list"): Settermethod for the commonBpsC2 slot.
- seqsC1<- signature(object = "Breakpoints", value = "list"): Setter-method
 for the seqsC1 slot.</pre>
- seqsC1 signature (object = "Breakpoints"): Getter-method for the seqsC1 slot.
- sqsC2<- signature(object = "Breakpoints", value = "list"): Setter-method
 for the seqsC2 slot.</pre>
- seqsC2 signature(object = "Breakpoints"): Getter-method for the seqsC2 slot.
- [signature(x = "Breakpoints", i = "ANY", j = "ANY"): Subsetting a Breakpoints object.
- **length** signature (x = "Breakpoints"): Returns the number of breakpoints stored.

- **names<-** signature (x = "Breakpoints", value = "ANY"): Set the names of the breakpoints.
- **names** signature (x = "Breakpoints"): Get the names of the breakpoints.
- plotChimericReads signature(brpData = "Breakpoints"): Plot the structural variant and the chimeric reads covering its breakpoints.

```
summary signature(object = "Breakpoints"): Create a data frame summaring in-
formation about all breakpoints.
```

table signature(... = "Breakpoints"): Create a frequency table of cluster sizes.

Author(s)

Hans-Ulrich Klein, Christoph Bartenhagen

See Also

filterChimericReads, detectBreakpoints, mergeBreakpoints, plotChimericReads

MapperSet-class

Class to Contain GS Reference Mapper Output

Description

Container to store data imported from a project of Roche's GS Reference Mapper Software. It stores all information into a Biobase ExpressionSet.

Objects from the Class

Objects can be created by calls of the form MapperSet (filename). While filename is a vector containing all sample directories (i.e. directories that contain the files "mapping/454HCDiffs.txt" and "mapping/454NewblerMetrics.txt").

Slots

- assayData: Object of class AssayData. Contains the number of reads with the respective difference and the total coverage for every variant in forward and reverse direction.
- featureData: Object of class AnnotatedDataFrame. Contains information about the type, location and reference of each variant. If available, it shows further Ensembl variant-ids for known SNPs.
- phenoData: Object of class AnnotatedDataFrame. By default, the phenoData contains the accession number of every sample.
- variantFilterPerc: Object of class numeric. Contains a threshold to display only those variants, whose coverage (in percent) in forward and reverse direction in at least one sample is higher than this filter value. See setVariantFilter for details about setting this value.
- variantFilter: Object of class character. Contains a vector of variant names whose coverage (in percent) in forward and reverse direction in at least one sample is higher than the filter value(s) in variantFilterPerc.
- dirs: Object of class character. Based on a directory given at instantiation of the object, it contains a vector of several directories containing all relevant GS Mapper project files.
- experimentData: Object of class MIAME. Contains details of the experiment.
- annotation: Object of class character Label associated with the annotation package used in the experiment.
- protocolData: Object of class AnnotatedDataFrame. Contains additional information about the samples.
- .__classVersion__: Object of class Versions. Remembers the R and R453Toolbox version numbers used to created the MapperSet instance.

MapperSet 9

Extends

Class eSet, directly. Class VersionedBiobase, by class "eSet", distance 2. Class Versioned, by class "eSet", distance 3.

Methods

setVariantFilter(object): Sets the filter to display only those variants, whose coverage (in percent) in forward and reverse direction in at least one sample is higher than the given value.

getVariantPercentages(object) Computes the coverage for every variant over all reads (forward and/or reverse) and for each sample.

annotateVariants(object): Annotates given genomic variants. See annotateVariants for details

htmlReport(object): Exports all (filtered) variant data into a html report. See htmlReport for details

getReadStatus(object): Reads the file "454ReadStatus.txt" in the GSM project directory which contains information about the alignment of each read (chr, pos, strand, etc.)and returns its content in a dataframe.

Author(s)

Christoph Bartenhagen

See Also

AVASet, annotateVariants, htmlReport, setVariantFilter, getVariantPercentages

Examples

```
# sum up class structure
showClass("MapperSet")

# load a GS Mapper dataset containing 3 samples and 111 variants
data(mapperSetExample)
mapperSetExample

# show contents of assay, feature and pheno data
assayData(mapperSetExample)
fData(mapperSetExample)
pData(mapperSetExample)
```

MapperSet

Creating a MapperSet

Description

This function imports a project of Roche's GS Reference Mapper Software. It stores all information into an instance of the Biobase ExpressionSet.

Usage

```
MapperSet (dirs, samplenames)
```

10 MapperSet

Arguments

dirs A character vector containing all sample directories (i.e. directories that contain

the files "mapping/454HCDiffs.txt" (required), "mapping/454ReadStatus.txt" (op-

tional), "mapping/454NewblerMetrics.txt"(optional)).

samplenames A character vector containing samplenames. The order and number of sample-

names must be consistent with the filenames to ensure the correctness of the MapperSet. If no samplenames are given, the filenames are used for naming.

Details

An instance of the MapperSet is derived from the Biobase eSet and thus structured into

1. assayData

variantForwCount/variantRevCount: Contain the number of reads with the respective difference in forward/reverse direction.

totalForwCount/totalRevCount: Contain the total coverage for every variant in forward/reverse direction.

2. featureData

chromosome, start/end: Give the location of each variant.

referenceBases/variantBase: Show the bases changed in each variant. **regName:** The name of the region (gene) where the variant is located.

knownSNP: Lists Ensembl variant-ids for known SNPs (if any).

3. phenoData

By default, the phenoData contains the accession number of every sample.

Value

An instance of the MapperSet.

Author(s)

Christoph Bartenhagen

See Also

AVASet-class

Examples

```
# load a GS Mapper dataset containing 3 samples and 111 variants
data(mapperSetExample)
mapperSetExample
```

SFFContainer-class 11

SFFContainer-class Class "SFFContainer"

Description

This class is a container for data from files in Roche's Standard Flowgram Format (SFF).

Objects from the Class

Objects can be created by calls of the form new ("SFFContainer", ...). Usually, objects will be created by calling the readSFF method on a file in SFF format.

Slots

- name: Object of class "character" containing the name of the file this SFFContainer was created from.
- flowgramFormat: Object of class "numeric" representing the format used to encode each of the flowgram values for each read. Currently, only one flowgram format has been adopted and is coded by the value 1.
- flowChars: Object of class "character" containing the array of nucleotide bases ('A', 'C', 'G', 'T') that correspond to the nucleotides used for each flow of each read.
- keySequence: Object of class "character" representing the nucleotide bases of the key sequence used for these reads.
- clipQualityLeft: Object of class "numeric" representing the position of the first base after the clipping point for an attached quality sequence for each read. If only a combined (quality+adapter) clipping position is computed it should be stored in clipQualityLeft. If no clipping value is computed the field is set to 0. The position values use 1-based indexing.
- clipQualityRight: Object of class "numeric" representing the position of the last base before the clipping point for an attached quality sequence for each read. If only a combined (quality+adapter) clipping position is computed it should be stored in clipQualityRight. If no clipping value is computed the field is set to 0. The position values use 1-based indexing.
- clipAdapterLeft: Object of class "numeric" representing the position of the first base after the clipping point for an attached adapter sequence for each read. If only a combined (quality+adapter) clipping position is computed it should be stored in clipQualityLeft. If no clipping value is computed the field is set to 0. The position values use 1-based indexing.
- clipAdapterRight: Object of class "numeric" representing the position of the last base before the clipping point for an attached adapter sequence for each read. If only a combined (quality+adapter) clipping position is computed it should be stored in clipQualityRight. If no clipping value is computed the field is set to 0. The position values use 1-based indexing.
- flowgrams: Object of class "list" containing the homopolymer stretch estimates for each flow using one list item for each read.
- flowIndexes: Object of class "list" containing the flow positions for each base in the called sequence, i.e. for each base, the position in the flowgram whose estimate resulted in that base being called. Each read has its own list item.
- reads: Object of class "QualityScaledDNAStringSet" containing the basecalled nucleotide sequences of each read together with the quality scores for each of the bases in the sequence using the standard -log10 probability scale.

12 SFFContainer-class

Methods

```
addRead signature(object = "SFFContainer", read = "SFFRead"): Adds an
  object of class SFFRead to the SFFContainer
```

- getRead signature(object = "SFFContainer", readname = "character"): Returns the read with the given name as an object of class SFFRead.
- clipAdapterLeft<- signature(object = "SFFContainer", value = "numeric"):
 Setter-method for the clipAdapterLeft slot.</pre>
- clipAdapterLeft signature(object = "SFFContainer"): Getter-method for the clipAdapterLeft slot.
- clipAdapterRight<- signature(object = "SFFContainer", value = "numeric"):
 Setter-method for the clipAdapterRight slot.</pre>
- clipAdapterRight signature(object = "SFFContainer"): Getter-method for the clipAdapterRight slot.
- clipQualityLeft<- signature(object = "SFFContainer", value = "numeric"):
 Setter-method for the clipQualityLeft slot.</pre>
- clipQualityLeft signature(object = "SFFContainer"): Getter-method for the clipQualityLeft slot.
- clipQualityRight<- signature(object = "SFFContainer", value = "numeric"):
 Setter-method for the clipQualityRight slot.</pre>
- clipQualityRight signature(object = "SFFContainer"): Getter-method for the clipQualityRight slot.
- name<- signature(object = "SFFContainer", value = "character"): Settermethod for the name slot.</pre>
- name signature(object = "SFFContainer"): Getter-method for the name slot.
- flowChars<- signature(object = "SFFContainer", value = "character"): Settermethod for the flowChars slot.</pre>
- flowgramFormat<- signature(object = "SFFContainer", value = "numeric"):
 Setter-method for the flowgramFormat slot.</pre>
- flowgramFormat signature(object = "SFFContainer"): Getter-method for the flowgramFormat slot.
- flowgrams<- signature(object = "SFFContainer", value = "list"): Setter-method
 for the flowgrams slot.</pre>
- flowIndexes<- signature(object = "SFFContainer", value = "list"): Settermethod for the flowIndexes slot.</pre>
- flowIndexes
 signature(object = "SFFContainer"): Getter-method for the flowIndexes
 slot.
- keySequence<- signature(object = "SFFContainer", value = "character"):
 Setter-method for the keySequence slot.</pre>
- keySequence signature(object = "SFFContainer"): Getter-method for the keySequence slot.
- reads<- signature(object = "SFFContainer", value = "QualityScaledDNAStringSet"):
 Setter-method for the reads slot.</pre>

SFFRead-class 13

```
reads signature(object = "SFFContainer"): Getter-method for the reads slot.
[ signature(x = "SFFContainer", i = "ANY", j = "ANY"): Subsetting a SFF-
Container object.
```

Author(s)

Christian Ruckert

See Also

```
readSFF, SFFRead
```

Examples

```
showClass("SFFContainer")
```

SFFRead-class

Class '"SFFRead"'

Description

This class is a container for a single read from files in Roche's Standard Flowgram Format (SFF).

Objects from the Class

Objects can be created by calls of the form new ("SFFRead", ...). Usually, objects will be created by calling the getRead method on an object of class SFFContainer.

Slots

name: Object of class "character" representing the name of the read.

read: Object of class "DNAString" containing the basecalled nucleotide sequence of the read.

- flowgramFormat: Object of class "numeric" representing the format used to encode each of the flowgram values for each read. Currently, only one flowgram format has been adopted and is coded by the value 1.
- flowChars: Object of class "character" containing the array of nucleotide bases ('A', 'C', 'G', 'T') that correspond to the nucleotides used for each flow of each read.
- keySequence: Object of class "character" representing the nucleotide bases of the key sequence used for these reads.
- clipQualityLeft: Object of class "numeric" representing the position of the first base after the clipping point for an attached quality sequence. If only a combined (quality+adapter) clipping position is computed it should be stored in clipQualityLeft. If no clipping value is computed the field is set to 0. The position values use 1-based indexing.
- clipQualityRight: Object of class "numeric" representing the position of the last base before the clipping point for an attached quality sequence. If only a combined (quality+adapter) clipping position is computed it should be stored in clipQualityRight. If no clipping value is computed the field is set to 0. The position values use 1-based indexing.

14 SFFRead-class

clipAdapterLeft: Object of class "numeric" representing the position of the first base after the clipping point for an attached adapter sequence. If only a combined (quality+adapter) clipping position is computed it should be stored in clipQualityLeft. If no clipping value is computed the field is set to 0. The position values use 1-based indexing.

- clipAdapterRight: Object of class "numeric" representing the position of the last base before the clipping point for an attached adapter sequence. If only a combined (quality+adapter) clipping position is computed it should be stored in clipQualityRight. If no clipping value is computed the field is set to 0. The position values use 1-based indexing.
- flowgram: Object of class "numeric" containing the homopolymer stretch estimates for each
- flowIndexes: Object of class "numeric" containing the flow positions for each base in the called sequence, i.e. for each base, the position in the flowgram whose estimate resulted in that base being called.
- quality: Object of class "BString" containing the quality scores for each of the bases in the sequence, where the values use the standard -log10 probability scale.

Methods

- read<- signature(object = "SFFRead", value = "DNAString"): Setter-method
 for the read slot.</pre>
- read signature(object = "SFFRead"): Getter-method for the read slot.
- flowChars<- signature(object = "SFFContainer", value = "character"): Settermethod for the flowChars slot.</pre>
- flowgramFormat<- signature(object = "SFFContainer", value = "numeric"):
 Setter-method for the flowgramFormat slot.</pre>
- flowgramFormat signature(object = "SFFContainer"): Getter-method for the flowgramFormat slot.
- keySequence<- signature(object = "SFFContainer", value = "character"):
 Setter-method for the keySequence slot.</pre>
- keySequence signature(object = "SFFContainer"): Getter-method for the keySequence slot.
- clipAdapterLeft<- signature(object = "SFFRead", value = "numeric"): Settermethod for the clipAdapterLeft slot.</pre>
- clipAdapterLeft signature(object = "SFFRead"): Getter-method for the clipAdapterLeft slot.
- clipAdapterRight<- signature(object = "SFFRead", value = "numeric"): Settermethod for the clipAdapterRight slot.</pre>
- clipAdapterRight signature(object = "SFFRead"): Getter-method for the clipAdapterRight slot.
- clipQualityLeft<- signature(object = "SFFRead", value = "numeric"): Settermethod for the clipQualityLeft slot.</pre>
- clipQualityRight<- signature(object = "SFFRead", value = "numeric"): Settermethod for the clipQualityRight slot.</pre>

alignShortReads 15

```
clipQualityRight signature(object = "SFFRead"): Getter-method for the clipQuali-
tyRight slot.
```

flowgram<- signature(object = "SFFRead", value = "numeric"): Setter-method
 for the flowgram slot.</pre>

flowgram signature(object = "SFFRead"): Getter-method for the flowgram slot.

flowIndexes<- signature(object = "SFFRead", value = "numeric"): Setter-method
 for the flowIndexes slot.</pre>

flowIndexes signature (object = "SFFRead"): Getter-method for the flowIndexes slot.

name<- signature(object = "SFFRead", value = "character"): Setter-method
 for the name slot.</pre>

name signature(object = "SFFRead"): Getter-method for the name slot.

quality<- signature(object = "SFFRead", value = "BString"): Setter-method
 for the quality slot.</pre>

quality signature (object = "SFFRead"): Getter-method for the quality slot.

Author(s)

Christian Ruckert

See Also

```
readSFF, SFFContainer
```

Examples

showClass("SFFRead")

alignShortReads

Exact alignment of DNA sequences against a reference

Description

This method aligns given sequences against a given reference genome using the matchPDict method. Only exact (no errors) and unique matches are returned.

Usage

alignShortReads(object, bsGenome, seqNames, ensemblNotation)

Arguments

object The reads that should be aligned agiven either as a DNAStringSet or a AVASet

instance. In the latter case the reference sequences are extracted and aligned.

bsGenome instance providing the reference sequences.

seqNames The names of the sequences in bsGenome that should be used. If omitted, all

reference sequences are used.

ensemblNotation

If set to TRUE, "chr" is removed from the reference sequences' names in the

returned alignment. Default value is FALSE.

16 annotate Variants

Details

All reads are aligned against the reference and its reverse complement. If the reads are not in 5' to 3' orientation, they should be reversed before. Note that only exact and unique alignments are reported. Use matchPDict directly for more flexibility.

Value

An object of class AlignedRead or a AVASet instance.

Author(s)

Hans-Ulrich Klein

See Also

```
matchPDict, DNAStringSet, AlignedRead, AVASet
```

Examples

```
library("BSgenome.Scerevisiae.UCSC.sacCer2")
reads = DNAStringSet(c(
   "CCGTTCAAAGAGCCCTTGGCCCATAATCCACCGGTT",
   "ATCCTGCCACAGGAGTCCATGGAGGTTTCGCCA"))
alignShortReads(reads, Scerevisiae, seqNames="chrIII")
```

annotateVariants

Adds genomic information to variants

Description

This method annotates given genomic variants (mutations). Annotation includes affected genes, exons and codons. Resulting amino acid changes are returned as well as dbSNP identifiers, if the mutation is already known. All information is fetched from Ensembl via biomaRt using the datasets hsapiens_gene_ensembl and hsapiens_snp.

Usage

```
annotateVariants(object, bsGenome)
```

Arguments

object A data frame storing variants or an instance of AVASet/MapperSet or a data

frame (see details).

bsGenome An object of class BSGenome giving the genome to be used as reference se-

quence to calculate amino acid changes. This argument is only applicable when object is of type MapperSet. Default is 'BSgenome.Hsapiens.UCSC.hg19'.

Note that the genome should fit to the Ensembl annotation.

Details

If a data frame is given, the following columns must be present:

assayDataAmp 17

start	genomic start position in the current Ensembl genome
end	genomic end position in the current Ensembl genome
chromosome	chromosome in ensembl notation (i.e. "1", "2",, "Y")
strand	"+" or "-" relative to the nucleotide bases given below
D C	

seqRef reference sequence

seqMut sequence of the observed variant

seqSur reference sequence extended for 3 bases in both directions

The rownames of the data frame are used as mutations' names (IDs). See examples for a properly defined data drame.

Value

An object of class AnnotatedVariants. Affected genes, transcripts and exon as well as known SNPs are stored in a list-like structure. See the documentation of class AnnotatedVariants—class for details.

Author(s)

Hans-Ulrich Klein

See Also

AnnotatedVariants-class, AVASet-class, MapperSet-class, htmlReport

Examples

```
variants = data.frame(
    start=c(106157528, 106154991,106156184),
    end=c(106157528, 106154994,106156185),
    chromosome=c("4", "4", "4"),
    strand=c("+", "+", "+"),
    seqRef=c("A", "ATAG", "---"),
    seqMut=c("G", "----", "ATA"),
    seqSur=c("TACAGAA", "TTTATAGATA", "AGC---TCC"),
    stringsAsFactors=FALSE)
rownames(variants) = c("snp", "del", "ins")
## Not run: annotateVariants(variants)
```

assayDataAmp

Access the amplicon data of an AVASet.

Description

Similar to assayData of the Biobase ExpressionSet, this function returns the assay data of the amplicon slot of an instance of the AVASet.

Usage

```
assayDataAmp(object)
```

18 avaSetExample

Arguments

```
object An link (AVASet-class) object.
```

Value

The assay data of the amplicon slot consists of a list of two data frames with the number of forward and reverse reads of all amplicons for each sample (see AVASet-class for details).

Author(s)

Christoph Bartenhagen

See Also

```
fDataAmp, featureDataAmp, AVASet-class
```

Examples

```
\# load an AVA dataset containing 6 samples, 4 amplicons and 259 variants data(avaSetExample)
```

```
# show contents of amplicon assay data
assayDataAmp(avaSetExample)
```

avaSetExample

Amplicon Variant Analyzer data import

Description

This is an example of an link {AVASet-class} object containing the output of Roche's Amplicon Variant Analyzer Software. It consists of 6 samples, 4 amplicons and 259 variants.

Usage

```
data(avaSetExample)
```

Format

Formal class 'AVASet'

Source

'Next-generation sequencing technology reveals a characteristic pattern of molecular mutations in 72.8 leukemia by detecting frequent alterations in TET2, CBL, RAS, and RUNX1' (Kohlmann A et al., J Clin Oncol. 2010 Aug 20;28(24):3858-65. Epub 2010 Jul 19)

Examples

```
data(avaSetExample)
avaSetExample
```

avaSetFiltered 19

avaSetFiltered

Amplicon Variant Analyzer data import

Description

This is an example of an link {AVASet-class} object containing the output of Roche's Amplicon Variant Analyzer Software. It consists of 6 samples, 4 amplicons and 4 variants. The variants were previously filtered according to the amplicon coverage (see setVariantFilter for details about filtering an AVASet object).

Usage

```
data(avaSetFiltered)
```

Format

Formal class 'AVASet'

Source

'Next-generation sequencing technology reveals a characteristic pattern of molecular mutations in 72.8 leukemia by detecting frequent alterations in TET2, CBL, RAS, and RUNX1' (Kohlmann A et al., J Clin Oncol. 2010 Aug 20;28(24):3858-65. Epub 2010 Jul 19)

Examples

data(avaSetFiltered)
avaSetFiltered

avaSetFiltered_annot

AVASet variant annotations

Description

These are example annotations for 4 variants of an AVASet (try data (avaSetFiltered) to retrieve the corresponding link{AVASet-class} object). The annotations include affected genes, exons and codons as well as resulting amino acid changes and dbSNP identifiers (if the mutation is already known).

Usage

```
data(avaSetFiltered_annot)
```

Format

Formal class 'Annotated Variants'

20 baseFrequency

Source

'Next-generation sequencing technology reveals a characteristic pattern of molecular mutations in 72.8 leukemia by detecting frequent alterations in TET2, CBL, RAS, and RUNX1' (Kohlmann A et al., J Clin Oncol. 2010 Aug 20;28(24):3858-65. Epub 2010 Jul 19)

Examples

data(avaSetFiltered_annot)

baseFrequency

Absolute And Relative Frequency Of The Four Bases.

Description

This function returns the absolute and the relative frequency of the four bases (A, C, G, T).

Usage

baseFrequency(object)

Arguments

object

An object of class DNAStringSet, ShortRead or SFFContainer.

Details

This function makes use of the alphabetFrequency function from package Biostrings.

Value

A data.frame with two columns containing the absolute and relative frequencies respectively and six rows, one for each of the four bases (A, C, G, T), one for other symbols contained in the reads and one summarizing the five aforementioned rows.

Author(s)

Christian Ruckert

baseQualityHist 21

baseQualityHist Plot A Histogram Of The Base Qualities.

Description

Create a histogram based on the quality of every single base from all sequences.

Usage

```
baseQualityHist(object, xlab="Quality score", ylab="Number of bases", col="fir
...)
```

Arguments

object An object of class QualityScaledDNAStringSet, ShortReadQ or SFFContainer.

xlab The X axis label.

ylab The Y axis label.

col The plotting color.

breaks The number of breaks in the histogram (see 'hist').

... Arguments to be passed to methods, such as graphical parameters (see 'par').

Author(s)

Christian Ruckert

baseQualityStats Statistics Of Base Quality

Description

This function returns mean, minimum, maximum and standard deviation of the base quality scores over all sequences.

Usage

```
baseQualityStats(object)
```

Arguments

object An object of class QualityScaledDNAStringSet, ShortReadQ or SFFContainer.

Value

A numeric vector with four slots: mean, min, max, sd.

Author(s)

Christian Ruckert

22 calculateTiTv

breakpoints

Putative breakpoints of chimeric reads

Description

This example holds two consensus (pathogenic and reciproce) breakpoints of 12 chimeric reads indicating an inversion on chromosome 16. The Breakpoints object gives access to the breakpoint locations as well as alignment information for each of the 12 reads.

Usage

```
data(breakpoints)
```

Format

Formal class 'Breakpoints'

Source

'Targeted next-generation sequencing detects point mutations, insertions, deletions, and balanced chromosomal rearrangements as well as identifies novel leukemia-specific fusion genes in a single procedure' (Leukemia, submitted)

Examples

```
data(breakpoints)
```

calculateTiTv

Calculate transition transversion ratio

Description

When many point mutations are detected, the ration of transitions to transversions can be used as quality measure to assess the number of false positives.

Usage

```
## S4 method for signature 'AVASet'
calculateTiTv(object)
  ## S4 method for signature 'MapperSet'
calculateTiTv(object)
```

Arguments

object

An instance of AVASet or MapperSet storing the detected variants.

Details

For more information about the Ti/Tv ratio see http://www.broadinstitute.org/gsa/wiki/index.php/QC_Methods

captureArray 23

Value

A list with two elements: A substitution matrix summarizing all observed substitutions and the transition/transversion ratio.

Author(s)

Hans-Ulrich Klein

Examples

```
data(avaSetExample)
ava = setVariantFilter(avaSetExample, c(0.03, 0.03))
calculateTiTv(ava)
```

captureArray

Custom capture array design

Description

Design of a custom Roche NimbleGen 385k capture array. The array captures short segments corresponding to all exon regions of 92 distinct target genes (genome build hg19). In addition, contiguous genomic regions were represented for three additional genes, i.e. CBFB, MLL, and RUNX1.

Usage

```
data(captureArray)
```

Format

Formal class 'CompressedIRangesList'

Source

'Targeted next-generation sequencing detects point mutations, insertions, deletions, and balanced chromosomal rearrangements as well as identifies novel leukemia-specific fusion genes in a single procedure' (Leukemia, submitted)

Examples

```
data(captureArray)
```

24 complexity.dust

complexity.dust	Sequence Complexity Using The DUST Algorithm

Description

This function evaluates the sequence complexity using the DUST algorithm.

Usage

```
complexity.dust(object, xlab="Complexity score (0=high, 100=low)", ylab="Number xlim=c(0, 100), col="firebrick1", breaks=100, ...)
```

Arguments

object	An object of class DNAStringSet, ShortRead or SFFContainer.
xlab	The X axis label.
ylab	The Y axis label.
xlim	The limits of the X axis.
col	The plotting color.
breaks	The number of breaks in the histogram (see 'hist').
	Arguments to be passed to methods, such as graphical parameters (see 'par').

Details

The complexity score is based on how often different trinucleotides occur and is scaled between 0 and 100. A sequence of homopolymer repeats (e.g. TTTTTTTTT) has a score of 100, of dinucleotide repeats (e.g. TATATATATA) has a score around 49, and of trinucleotide repeats (e.g. TAGTAGTAG) has a score around 32. Scores above seven can be considered low-complexity.

Value

A numeric vector containing the complexity score for each sequence.

Author(s)

Christian Ruckert

References

Schmieder R. (2011) Quality control and preprocessing of metagenomic datasets. *Bioinformatics*, 2011 Mar 15;27(6):863-4.

complexity.entropy 25

complexity.entropy Sequence Complexity Using The Shannon-Wiener Algorithm

Description

This function evaluates the sequence complexity using the Shannon-Wiener Algorithm.

Usage

```
complexity.entropy(object, xlab="Complexity score (0=low, 100=high)", ylab="Nu
xlim=c(0, 100), col="firebrick1", breaks=100, ...)
```

Arguments

object	An object of class DNAStringSet, ShortRead or SFFContainer.
xlab	The X axis label.
ylab	The Y axis label.
xlim	The limits of the X axis.
col	The plotting color.
breaks	The number of breaks in the histogram (see 'hist').
	Arguments to be passed to methods, such as graphical parameters (see 'par').

Details

The entropy approach evaluates the entropy of trinucleotides in a sequence. The entropy values are scaled from 0 to 100 and lower entropy values imply lower complexity. A sequence of homopolymer repeats (e.g. TTTTTTTTTT) has an entropy value of 0, of dinucleotide repeats (e.g. TATATATATA) has an entropy value around 16, and of trinucleotide repeats (e.g. TAGTAGTAG) has an entropy value around 26. Scores below 70 can be considered low-complexity.

Value

A numeric vector containing the complexity score for each sequence.

Author(s)

Christian Ruckert

References

Schmieder R. (2011) Quality control and preprocessing of metagenomic datasets. *Bioinformatics*, 2011 Mar 15;27(6):863-4.

26 coverageOnTarget

convertCigar

Basic functions for CIGAR strings

Description

These are temporary methods, that are likely to be replaced by methods from the Rsamtools package in near future.

Usage

```
extendedCIGARToList(cigars)
listToExtendedCIGAR(cigarList)
```

Arguments

cigars A character vector with CIGAR strings.

cigarList A list of converted CIGAR strings as produced by extendedCIGARToList

coverageOnTarget

Computes the coverage restricted to the target region.

Description

This method computes the approximate coverage of each base in a given region.

Usage

```
coverageOnTarget(alnReads, targetRegion)
```

Arguments

alnReads A list as returned by scanBam storing aligned reads.

targetRegion The target region as a RangesList. The chromosome names must fit to the chromosome names used in the alignment information of the given reads.

Details

The detailed alignment information given by the CIGAR strings in .bam files are ignored by the function. Instead, it is assumed that the whole read alignes to the reference without indels. This is often not true for longer read (e.g. generated with Roche 454 Sequencing), but saves computation time.

Value

A list of the same length as the alnReads argument. Each list element is an integer vector of the same length as the target region (in bases) and stores the coverage generated by the reads from the corresponding list element of alnReads.

demultiplexReads 27

Author(s)

Hans-Ulrich Klein

See Also

scanBam

Examples

```
library(Rsamtools)
bamFile = system.file("extdata", "StructuralVariantDetection", "bam", "N01.bam", package=
bam = scanBam(bamFile)
region = RangesList("11"=IRanges(start=118307205, end=118395936))
cov = coverageOnTarget(bam, region)
```

demultiplexReads

Performs MID/Multiplex filtering

Description

Roche's Genome Sequencer allows to load two or more samples on one region. To allocate sequences to samples, each sample has a unique multiplex sequence. The multiplex sequence should be the prefix of all sequences from that sample. This method demultiplexes a given set of sequences according to the given multiplex sequences (MIDs).

Usage

```
## S4 method for signature 'XStringSet, XStringSet, numeric, logical'
demultiplexReads(reads, mids, numMismatches, trim)
```

Arguments

reads A DNAStringSet instance that contains reads starting with MIDs

mids A DNAStringSet instance that contains the MIDs

numMismatches

The maximal number of mismatches allowed, default 2.

trim Whether the MIDs should be cutted-out, default TRUE

Details

All given MIDs must have the same length. The algorithm computes the number of mismachtes for each MID. The read is assigned to the MID with the lowest number of mismaches. If two or more MIDs have the same number of mismachtes, or if the number of mismachtes is greater than the given argument numMismachtes, the read is not assigned to any MID. The default number of allowed mismatches is 2.

Value

demultiplexReads returns a list with one DNAStringSet instance for each MID.

28 detectBreakpoints

Author(s)

Hans-Ulrich Klein

See Also

```
genomeSequencerMIDs, DNAStringSet
```

Examples

```
library(Biostrings)
  mids = genomeSequencerMIDs(c("MID1", "MID2", "MID3"))
  reads = DNAStringSet(c(
      paste(as.character(mids[["MID1"]]), "A", sep=""),
      paste(as.character(mids[["MID1"]]), "AAA", sep=""),
      paste(as.character(mids[["MID2"]]), "AAA", sep="")))
  demultiplexReads(reads, mids)
```

detectBreakpoints Clustering and consensus breakpoint detection for chimeric reads

Description

Given a set of chimeric reads, this methods computes all putative breakpoints. First, chimeric reads are clustered such that all reads spanning the same breakpoint form a cluster. Then, a consensus breakpoint sequence and breakpoint position is computed for each cluster.

Usage

```
detectBreakpoints(chimericReads, bpDist=100, minClusterSize=4, removeSoftClips=T
```

Arguments

chimericReads

A list storing chimeric reads as returned by filterChimericReads. The list must have the format as defined by the scanBam method.

bpDist

The maximum distance in base pairs between the breakpoints of two chimeric reads at which the reads are merge to a cluster.

minClusterSize

Cluster whose size is below minClusterSize are be excluded from breakpoint detection.

removeSoftClips

If true, soft-clipped bases at the beginning or the end of a sequence are removed (see details below).

bsGenome

A bsGenome instance providing the reference sequences. If missing, the library BSgenome. Hsapiens. UCSC.hg19 is used by default.

detectBreakpoints 29

Details

This method is usually invoked after calling filterChimericReads and before calling mergeBreakpoints. It first forms clusters of chimeric reads (reads with exactly two local alignments) that span the same breakpoint and than computes a consensus breakpoint sequence for each cluster.

To carry out a hierarchical clustering, a measure for the distance between two chimeric reads must be defined. If reads span different chromosomes, their distance is set to infinity. The strand information of the local alignments may also indicate that two chimeric reads do not span the same breakpoint even if they span the same chromosomes. For example, the first reads has two local alignments on the positive strand whereas the second read has one local alignment on the positive strand and the other on the negative strand. In this case, the distance is set to inifinty, too. Finally, the distance measure distinguishes between the two breakpoints (sometimes called the pathogenic and the reciproce breakpoint) that originate from the same structual variant. The distance between a read from the pathogenic and a read from the reciproce breakpoint is infinity so that two different clusters will emerge. These two related breakpoints can be merge later using the mergeBreakpoints method. We observed that the breakpoints of these two cases often differ by a few ten or even a few hundred basepairs.

If the chromosome and strand information between two reads x and y are coherent, the Euclidian distance is used:

$$d(x,y) = (bp(x,ChrA) - bp(y,ChrA))^2 + (bp(x,ChrB) - bp(y,ChrB))^2$$

where bp gives the coordinates of the breakpoint for the given read and chromosome. Hierarchical clustering is applied with complete linkage and the dendrogram is cutted at a height of bpDist to obtain the final clusters. The bpDist argument does usually not influence the result, because we observed that reads spanning the same breakpoint have very little variation (only a few base pairs) in their local alignments due to sequencing errors or due to ambiguity caused by same/similar sequence of both chromosome near the breakpoint.

Although the given set of reads may belong to the same chimeric DNA, their individual breakpoints may differ in a few base pairs. Furthermore, a single read may have more than one possible breakpoint if a (small) part of the read was aligned to both parts.

The following step determines a consensus breakpoint for each cluster. It uses the supplied bsGenome to construct a chimeric reference sequence for all possible breakpoints over all reads within each cluster. After the reads were realigned to the chimeric reference sequences, the one that yields the highest alignment score is taken to represent best the chimeric DNA and its breakpoints.

As a preprocessing step, detectBreakpoints offers to remove soft clips occuring after the alignment:

Some reads may contain soft-clipped bases (e.g. linker sequences) at the beginning of the first part of the read or at the end of the second part. By default, detectBreakpoints removes these unaligned subsequences and adjusts the cigar string, the sequence, the sequence width (qwidth) and the local start/end coordinates.

Value

detectBreakpoints returns an object of class breakpoints, which is a list of breakpoint clusters, which gives access to all alignments and consensus breakpoints:

seqs

This IRanges DataFrame is mainly a rearranged version of the alignment input in chimericReads. In addition, it shows the corresponding breakpoints and local alignment coordinates.

30 detectBreakpoints

A dataframe listing the breakpoints for both parts of the chimeric reference, the associated chromosome, strand and the reference sequence itself, including positions "localStart"/"localEnd" indicating which part of the reference belongs to which breakpoint.

CommonAlign An object of class PairwiseAlignedFixedSubject of the Biostrings package that contains the alignment to the (best) consensus reference sequence.

alignedReads On the basis of commonAlign and commonBps, alignedReads is an instance of class AlignedRead containing all aligned reads including their associated chromosomes, strands, and positions. Since the reference is a chimeric

sequence each read has two chromosome and two strand entries.

Author(s)

Hans-Ulrich Klein, Christoph Bartenhagen

See Also

filterChimericReads mergeBreakpoints plotChimericReads

Examples

```
# Construct a small example with three chimeric reads
# (=6 local alignments) in bam format as given by
# aligners such as BWA-SW.
# The first two reads originate from the same case but
# from different strands. The third read originate from
# the reciprocal breakpoint.
library("BSgenome.Scerevisiae.UCSC.sacCer2")
bamReads = list()
bamReads[[1]] = list(
    qname=c("seq1", "seq1", "seq2", "seq2", "seq3", "seq3"),
    flag = as.integer(c(0, 0, 16, 16, 0, 0)),
    rname = factor(c("II", "III", "III", "II", "III", "II")),
    strand = factor(c("+", "+", "-", "-", "+", "+")),
    pos = as.integer(c(99951, 200000, 200000, 99951, 199950, 100001)),
    qwidth = as.integer(c(100, 100, 100, 100, 100, 100)),
    cigar = c("50M50S", "50S50M", "50S50M", "50M50S", "50M50S", "50S50M"),
    seq = DNAStringSet(c(
        paste(substr(Scerevisiae$chrII, start=99951, stop=100000),
              substr(Scerevisiae$chrIII, start=200000, stop=200049),
              sep=""),
        paste(substr(Scerevisiae$chrII, start=99951, stop=100000),
              substr(Scerevisiae$chrIII, start=200000, stop=200049),
              sep=""),
        paste(substr(Scerevisiae$chrIII, start=200000, stop=200049),
              substr(Scerevisiae$chrII, start=99951, stop=100000),
              sep=""),
        paste(substr(Scerevisiae$chrIII, start=200000, stop=200049),
              substr(Scerevisiae$chrII, start=99951, stop=100000),
              sep=""),
        paste(substr(Scerevisiae$chrIII, start=199950, stop=199999),
              substr(Scerevisiae$chrII, start=100001, stop=100050),
              sep=""),
        paste(substr(Scerevisiae$chrIII, start=199950, stop=199999),
              substr(Scerevisiae$chrII, start=100001, stop=100050),
```

dinucleotideOddsRatio 31

```
sep="")))
)

bps = detectBreakpoints(bamReads, minClusterSize=1, bsGenome=Scerevisiae)
summary(bps)
table(bps)
mergeBreakpoints(bps)
```

dinucleotideOddsRatio

Dinucleotide Odds Ratio

Description

This function calculates the dinucleotide odds ratio for each of the sixtheen possible dinucleotides.

Usage

```
dinucleotideOddsRatio(object, xlab="Under-/over-representation of dinucleotide
  col="firebrick1", ...)
```

Arguments

object An object of class DNAStringSet, ShortRead or SFFContainer.

xlab The X axis label.

col The plotting color.

Arguments to be passed to methods, such as graphical parameters (see 'par').

Details

The dinucleotide odds ratio assigns a value between 0 and 2 to each of the sixtheen possible dinucleotides (AA, AC, AG, AT, CA, CC, CG, CT, GA, GC, GG, GT, TA, TC, TG, TT). For values below 1 the dinucleotide is under-represented compared to the randomly expected frequency of this dinucleotide in a sequence of the given length and with the given frequencies of the four nucleotides (A, C, G, T). For values above 1 this dinucleotide is over-represented.

Value

A matrix with sixtheen columns, one for each dinucleotide, containing the dinucleotide odds ratio values for each sequence in a seperate row.

Author(s)

Christian Ruckert

References

Schmieder R. (2011) Quality control and preprocessing of metagenomic datasets. *Bioinformatics*, 2011 Mar 15;27(6):863-4.

32 featureDataAmp

fDataAmp

Access the amplicon data of an AVASet.

Description

Similar to fData of the Biobase ExpressionSet, this function returns the feature data of the amplicon slot of an instance of the AVASet.

Usage

```
fDataAmp(object)
```

Arguments

object

An link{AVASet-class} object.

Value

The feature data of the amplicon slot contains the names, primers, start/end positions and reference sequences of all amplicons (seeAVASet-class for details). It returns a data frame.

Author(s)

Christoph Bartenhagen

See Also

featureDataAmp, assayDataAmp, AVASet-class

Examples

```
# load an AVA dataset containing 6 samples, 4 amplicons and 259 variants
data(avaSetExample)
avaSetExample

# show contents amplicon feature data
fDataAmp(avaSetExample)
```

featureDataAmp

Access the amplicon data of an AVASet

Description

Similar to featureData of the Biobase ExpressionSet, this function returns the feature data and feature meta of the amplicon slot of an instance of the AVASet.

Usage

```
featureDataAmp(object)
```

filterChimericReads 33

Arguments

object An link{AVASet-class} object.

Value

The feature data of the amplicon slot contains the names, primers, start/end positions and reference sequences of all amplicons (see AVASet-class for details). The returned object is of class AnnotatedDataFrame.

Author(s)

Christoph Bartenhagen

See Also

fDataAmp, assayDataAmp, AVASet-class,

Examples

```
\# load an AVA dataset containing 6 samples, 4 amplicons and 259 variants data(avaSetExample) avaSetExample
```

show contents amplicon feature data
featureDataAmp(avaSetExample)

filterChimericReads

Extract chimeric reads and apply filtering steps to remove artificial chimeric reads.

Description

Chimeric reads may be caused by sequencing a chromosomal aberration or by technical issues during sample preparation. This method implements several filter steps to remove false chimeric reads.

Usage

```
## S4 method for signature 'list,RangesList,DNAString,numeric,numeric'
filterChimericReads(alnReads, targetRegion, linkerSeq, minDist, dupReadDist)
```

Arguments

alnReads	A list storing the aligned reads as produced by the function scanBam.
targetRegion	A object of class rangesList containing the target region of e.g. a used capture array. The parameter may be omitted in case of a non targeted sequencing approach.
linkerSeq	A linker sequence that was used during sample preparation. It may be omitted.
minDist	The minimum distance between two local alignments (see details), default 1000
dupReadDist	The maximum distance between the 5 prime start position of two duplicated reads (see details), default 1.

34 filterChimericReads

Details

The following filter steps are performed:

1. All chimeric reads with exactly two local alignments are extracted. Reads with more than two local alignments are discarded.

- 2. If the targetRegion argument is given, chimeric reads must have one local alignment at least overlapping the the target region. If both local alignments are outside the target region, the read is discarded.
- 3. If the linkerSeq argument is given, all chimeric reads that have the linker sequence between their local alignments are removed. When searching the linker sequence, 4 mismatches or indels are allowed and the linker sequence must not start or end within the first or last ten bases of the read. The function searches for the linkerSeq and for it's reverse complement.
- 4. Two local alignment of a read must have minDist reads between the alignments (if both alignment are on the same chromosome). Otherwise, the read seems to span a deletion and not a chromosomal aberration and is discarded.
- 5. Duplicated reads are removed. Two reads are duplicated, if the lie on the same strand and have the same 5 prime start position. Due to sequencing and alignment errors, the start position may vary for a maximum of dupReadDist bases. In case of duplicated reads, only the longest read is kept.

Reads passing all filtering steps are returned in the list structure as given by the alnReads argument (as derived from the scanBam method). A data frame with information about the number of reads that passed each filter is added to the list.

Value

A list containing only filtered chimeric reads. The list has the same structure like the given argument alnReads. Additionally, one element "log" with logging information of each filtering step is added.

Author(s)

Hans-Ulrich Klein

See Also

detectBreakpoints, mergeBreakpoints, Breakpoints-class, scanBam, sequenceCaptureLinke

Examples

```
library(Rsamtools)
bamFile = system.file("extdata", "StructuralVariantDetection", "bam", "N01.bam", package=
bam = scanBam(bamFile)
data(captureArray)
linker = sequenceCaptureLinkers("gSel3")[[1]]
filterReads = filterChimericReads(bam, targetRegion=captureArray, linkerSeg=linker)
```

flowgramBarplot 35

flowgramBarplot

Create A Barplot Of The Flow Intensities

Description

This function creates a barplot of the flow intensities with one bar for each nucleotide in the flow. With the height giving the measured intensity.

Usage

```
flowgramBarplot(x, range=c(0, length(flowgram(x))), xlab="Flow sequence",
   ylab="Flow intensity", col=c(A="black", C="red", G="blue", T="green"), ...)
```

Arguments

X	An object of class SFFRead.
range	Two positions between which the flows should be plotted.
xlab	The X axis label.
ylab	The Y axis label.
col	The colors in which the four nucleotids should be plotted.
	Arguments to be passed to methods, such as graphical parameters (see 'par').

Author(s)

Christian Ruckert

gcContent

Calculate The Overall GC-Content

Description

This function calculates the GC-content summarized over all sequences.

Usage

```
gcContent (object)
```

Arguments

object An object of class DNAStringSet, ShortRead or SFFContainer.

Details

```
The GC-content is calculated as follows: (\#G + \#C / \#G + \#C + \#A + \#T) * 100
Where \#G is the number of base G in all sequences.
```

36 gcPerPosition

Value

A numeric vector of length one containing the overall GC-content.

Author(s)

Christian Ruckert

gcContentHist GC-Content Histogram

Description

This function creates a histogram of the relative GC-content per sequence.

Usage

```
gcContentHist(object, xlab="GC content", ylab="Number of sequences", col="fire
breaks=50, ...)
```

Arguments

object An object of class DNAStringSet, ShortRead or SFFContainer.

xlab The X axis label.

ylab The Y axis label.

col The plotting color.

breaks The number of breaks in the histogram (see 'hist').

Arguments to be passed to methods, such as graphical parameters (see 'par').

Author(s)

Christian Ruckert

gcPerPosition GC-Content Per Position

Description

This function plots the GC-content frequency per base position.

Usage

```
gcPerPosition(object, range=0.95, type="1", col="blue", xlab="Position", ylab=
```

Arguments

object	An object of class DNAStringSet, ShortRead or SFFContainer.
range	An integer vector of length one or two. If length one only bases up to the percent quantile of read lengths defined by the given value are shown. A vector of length two gives the start and end base between which the GC-content is plotted.
type	The type of the plot (see 'plot').
col	The plotting color.
xlab	The X axis label.
ylab	The Y axis label.
	Arguments to be passed to methods, such as graphical parameters (see 'par').

Author(s)

Christian Ruckert

```
genomeSequencerMIDs
```

Retrieve GS multiplex sequences

Description

This method returns the standard multiplex sequences used by the Genome Sequence MID library kits.

Usage

Arguments

mid Character vector with multiplex sequences' IDs (MIDs)

Details

If the argument mid is omitted, all 14 available multiplex sequences are returned.

Value

genomeSequencerMIDs returns a DNAStringSet with the requested multiplex sequences.

Author(s)

Hans-Ulrich Klein

See Also

```
demultiplexReads
```

38 getAlignedReads

Examples

```
genomeSequencerMIDs()
genomeSequencerMIDs(c("MID1", "MID3"))
```

getAlignedReads

Import reads from an Amplicon Variant Analyzer project

Description

For a given AVASet, this function imports all aligned reads belonging to all (or some selected) amplicons of all samples.

Usage

```
getAlignedReads(object, amplicons, dir)
```

Arguments

object An instance of the link {AVASet-class}.

amplicons An (optional) character vector of amplicon names as mentioned in the amplicon

feature data (see fDataAmp).

dir Usually, the method tries to retrieve the path to the AVA project from the given

link (AVASet) object. However, if it fails to find the directory, dir can be

used to set the root directory of the AVA project.

Details

This function reports all reads for all samples together. If you want to get the reads for some samples individually, try subsetting your AVASet as in the examples below.

Value

One DNAStringSet that contains all aligned reads for all samples (eventually restricted to some given amplicons).

Author(s)

Christoph Bartenhagen

See Also

```
AVASet-class, fDataAmp
```

```
# load an AVA dataset containing 6 samples, 4 amplicons and 259 variants
data(avaSetExample)

# import all reads for amplicon "TET2_E11.04" of the first sample
avaProjectDir = system.file("extdata", "AVASet", package = "R453Plus1Toolbox")
alnReads = getAlignedReads(avaSetExample[, 1], dir=avaProjectDir, amplicons="TET2_E11.04"
show(alnReads)
```

getAminoAbbr 39

getAminoAbbr

Get amino acid abbreviations

Description

This function returns a table with amino acid names as first column and the according abbreviations as row names.

Usage

```
getAminoAbbr()
```

getVariantPercentages

Variant coverage

Description

This function computes the coverage for each variant (in forward and/or reverse direction) for all samples. The coverage is defined as the percentual amount of reads that cover a variant.

Usage

```
getVariantPercentages(object, direction="both")
```

Arguments

object An instance of class AVASet-class or MapperSet-class.

direction A character indicating the direction ("forward", "reverse" or "both").

Details

If the direction was set to "both", the percentages are computed over the sum of both directions. Otherwise it is computed only over the occurences in one direction (forward or reverse). The occurences can be accesses via assayData.

Value

 ${\tt getVariantPercentages}\ {\tt returns}\ {\tt a}\ {\tt data}\ {\tt frame}\ {\tt with}\ {\tt all}\ {\tt percentages/frequencies}\ {\tt for}\ {\tt all}\ {\tt samples}.$

Author(s)

Christoph Bartenhagen

See Also

```
setVariantFilter.
```

40 homopolymerHist

Examples

```
# load a (filtered) AVA dataset containing 6 samples, 4 amplicons and 4 variants
data(avaSetFiltered)
avaSetFiltered

# both directions
getVariantPercentages(avaSetFiltered, direction="both")
# this is equivalent to
(assayData(avaSetFiltered)[[1]] + assayData(avaSetFiltered)[[3]]) / (assayData(avaSetFiltered)
# forward direction only
getVariantPercentages(avaSetFiltered, direction="forward")
# this is equivalent to
assayData(avaSetFiltered)[[1]] / assayData(avaSetFiltered)[[2]]
# reverse direction only
getVariantPercentages(avaSetFiltered, direction="reverse")
# this is equivalent to
assayData(avaSetFiltered)[[3]] / assayData(avaSetFiltered)[[4]]
```

homopolymerHist

Create A Histogram Of The Homopolymer Stretches

Description

This function creates a histogram for the different lengths of the homopolymer stretches with one bar for each nucleotide in the flow. With the height giving the number of this homopolymer stretch in the flowgram.

Usage

```
\label{lower} homopolymerHist(x, range=c(0, length(flowgram(x))), xlab="Homopolymer length", ylab="Number of homopolymers", col=c(A="black", C="red", G="blue", T="green", col=c(A="black", C="red", G="blue", C="red", C="re
```

Arguments

2	X	An object of class SFFRead.
	range	Two positions between which the flows should be plotted.
2	ĸlab	The X axis label.
7	ylab	The Y axis label.
(col	The colors in which the four nucleotids should be plotted.
		Arguments to be passed to methods, such as graphical parameters (see 'par').

Author(s)

Christian Ruckert

htmlReport 41

htmlReport	HTML-Report Builder for the AVASet and MapperSet	

Description

This function creates a HTML variant and quality report for a given AVASet or MapperSet instance.

Usage

```
htmlReport(object, annot, blocks=c(), transcripts=c(), sampleCols, minMut=3, dir
```

Arguments

object	An AVASet-class or MapperSet-class instance.
annot	An instance of class Annotated Variants you get by calling annotate Variants. If no such argument is supplied the data will be read from the Ensembl database automatically for all variants.
blocks	Character vector of block names for each variant. The variants will then be structured into several blocks. If no such list is supplied, the report consists of just one (big) table for all variants.
transcripts	Character vector containing Ensembl transcript-IDs that order the according entries on the transcript pages. Transcripts given in this argument will appear on top of the transcript page.
sampleCols	Character vector of column names of the sample data (phenoData) of the AVASet/MapperSet object to filter the sample output on the transcript pages. All columns will be listed if no such argument is given.
minMut	If the value of minMut is greater than zero, the report lists only variants, whose coverage for at least one sample is higher than minMut (percentage between 0 and 100).
dir	Character with the desired output directory. By defaultm the directory "HTML-Report" will be created in the current directory.
title	Heading for the first page with the variant information.

Details

The report is structured into two (MapperSet) or three (AVASet) parts containing variant and quality information:

- The main page sums up given variant information like the name, type, reference gene, position (see fData, annotateVariants).
 - Using the argument blocks, the main page can be individually structured by assigning a block name to each variant.
 - The main page can be further structured by samples. For a given AVASet object, every sample links to another short quality report showing only the amplicon coverage for this sample.
- Every variant on the main page links to a page with further details about the affected genes and transcripts (e.g. Ensembl gene-IDs, transcript-IDs, codon sequences, changes of amino acids (if coding)).
- Only in case of AVASet object: A quality report shows the coverage of every amplicon in forward and/or reverse direction. Further plots display the coverage by MID and PTP (if this information is given in the pheno data of the object).

42 mapperSetExample

Author(s)

Christoph Bartenhagen, Hans-Ulrich Klein, Christian Ruckert

See Also

```
annotateVariants.
```

Examples

```
# note: all examples save the report to the directory "htmlReportExample" in your current
# load a filtered AVA dataset containing 6 samples, 4 amplicons and 4 variants
# and its variant annotations
data("avaSetFiltered")
data("avaSetFiltered annot")
# create a full report showing all (unfiltered) information
htmlReport(avaSetFiltered, avaSetFiltered_annot, dir="htmlReportExample", title="htmlReport
# create a report that emphasizes on samples with variants covered by at least 50% of the
htmlReport(avaSetFiltered, avaSetFiltered_annot, dir="htmlReportExample", title="htmlReportExample",
# create a report that is structured by the reference genes
library("ShortRead")
refs = sapply(fData(avaSetFiltered)$referenceSeq, function(x)
    subset (pData(alignData(referenceSequences(avaSetFiltered))), pData(alignData(referenceSequences))
htmlReport(avaSetFiltered, avaSetFiltered_annot, dir="htmlReportExample", title="htmlReport
# create a report whose sample information only lists the sample ids
pData(avaSetFiltered)
sampleCols = "SampleID"
htmlReport(avaSetFiltered, avaSetFiltered_annot, dir="htmlReportExample", title="htmlReport
```

Description

This is an example of an link {MapperSet-class} object containing the output of Roche's GS Reference Mapper Software. It consists of 3 samples and 111 variants.

Usage

```
data(mapperSetExample)
```

Format

Formal class 'MapperSet'

Source

'Targeted next-generation sequencing detects point mutations, insertions, deletions, and balanced chromosomal rearrangements as well as identifies novel leukemia-specific fusion genes in a single procedure' (Leukemia, submitted)

mergeBreakpoints 43

Examples

data(mapperSetExample)
mapperSetExample

mergeBreakpoints

Identify and merge related breakpoints caused by the same variant.

Description

Structural variation like transversions or inversion cause two breakpoints. In the context of fusion genes, these are called the pathogenic breakpoint and the reciproce breakpoint. The method detectBreakpoints processes each breakpoint individually and does explicitly not put reads from the pathogenic and reciproce breakpoint into the same cluster. Hence, it is usually sensible to call this methods afterward to search for related pairs of breakpoints to gain more confidence about the existence of a structural variation.

Usage

mergeBreakpoints (breakpoints, maxDist, mergeBPs)

Arguments

breakpoints An object of class Breakpoints storing the breakpoints that will (potentially)

be merged.

maxDist The maximal distance in basepairs at which two breakpoints will be merged.

Default value is 1000.

mergeBPs An optional list of vectors of length two giving the breakpoints that should be

merged. If this argument is given, the method will not search for related break-

points.

Details

If the maxDist argument is given, the method compares each pair of breakpoints and checks, whether the two breakpoints may belong to the same structural variation. In addition to the spanned chromosomes, the orientation and the strand information of the reads spanning the breakpoints are also compared for this purpose. If chrosmosome, orientation and strand information of two breakpoints go well together, they will be merged, if the absolute distance of the breakpoints on chromosome A plus the absolute distance on chromosome B is smaller or equal to maxDist. If one breakpoint has more than one potential mate breakpoint for merging, it will be merged with the first candidate and a warning message is printed. The default value of maxDist is 1000.\

If the mergeBPs argument is given, the method will not search for related breakpoints but simply merge the given breakpoints. mergeBPs must be a list with vectors of length two that either contain the names of the indices of the breakpoints that should be merged.\

The arguments maxDist and mergeBPs cannot be given together. The given Breakpoints object must not contain breakpoints that have been merged before.

Value

An object of class Breakpoints storing merged and unmerged breakpoints.

44 nucleotideCharts

Author(s)

Hans-Ulrich Klein

See Also

```
detectBreakpoints, Breakpoints-class, plotChimericReads
```

Examples

```
# Load bam file and filter chimeric reads
library(Rsamtools)
bamFile = system.file("extdata", "StructuralVariantDetection", "bam", "N01.bam", package=bam = scanBam(bamFile)
data(captureArray)
linker = sequenceCaptureLinkers("gSel3")[[1]]
filterReads = filterChimericReads(bam, targetRegion=captureArray, linkerSeq=linker)

# detect breakpoints of size >= 3
breakpoints = detectBreakpoints(filterReads, minClusterSize=3)
table(breakpoints)
summary(breakpoints)

# merge breakpoints
breakpoints = mergeBreakpoints(breakpoints)
summary(breakpoints)
```

nucleotideCharts Nucleotide Charts

Description

This function plots the relative frequency of the four bases for each position in the sequences.

Usage

```
nucleotideCharts(object, range=0.95, linetypes=c(A="l", C="l", G="l", T="l", N
    linecols=c(A="black", C="red", G="blue", T="green", N="grey70"), xlab="Posit
    ylab="Frequency", ...)
```

Arguments

object	An object of class DNAStringSet, ShortRead or SFFContainer.
range	An integer vector of length one or two. If length one only bases up to the percent quantile of read lengths defined by the given value are shown. A vector of length two gives the start and end base between which the nucleotide frequencies are plotted.
linetypes	The line types used for the four nucleotids + N, see (see 'par')
linecols	The colors in which the four nucleotids + N should be plotted.
xlab	The X axis label.
ylab	The Y axis label.
	Arguments to be passed to methods, such as graphical parameters (see 'par').

plotAmpliconCoverage 45

Author(s)

Christian Ruckert

```
plotAmpliconCoverage
```

Creates a plot visualizing the number of reads per amplicon

Description

A function for visualizing the number of reads per amplicon or per MID / pico titer plate.

Usage

```
## S4 method for signature 'AVASet, character, logical'
plotAmpliconCoverage(avaSet, type="amplicon", bothDirections=TRUE, cex.names=0.8
```

Arguments

avaSet An instance of AVASet.

type A character vector specifying the type of plot.

bothDirections

A logical value determining whether the plot sums forward and reverse reads or

shows them separately.

cex.names Font size of the amplicon name labels.

cex.axis Font size of axes' labels.

las Orientation of amplicon name labels.

col Colors used in the plot.

... Arguments to be passed to methods, such as graphical parameters (see 'par').

Details

If the argumnet type is "amplicon", the number of reads for each amplicon are visualized. In case of a AVASet with one sample, a barplot with one bar for each amplicon is created. In case of more than one sample, a boxplot with one box for each amplicon is plotted. If type is "mid", a boxplot with one box for each MID is created. If type is "ptp", a boxplot with one box for each pico titer plate is created.

Author(s)

Hans-Ulrich Klein

See Also

AVASet

```
## Not run: data(avaSetExample)
plotAmpliconCoverage(avaSetExample)
plotAmpliconCoverage(avaSetExample[,1])
## End(Not run)
```

46 plotChimericReads

```
plotChimericReads Plots chimeric reads
```

Description

This function plots a given set of aligned chimeric reads along a reference sequence. It plots the breakpoints of translations or inversions and marks deletions, insertions and mismatches. Optionally, it displays all base pairs in a given region around the breakpoint.

Usage

Arguments

brpData	A Breakpoints object containing the consensus breakpoint of all reads and the consensus reference sequence as returned by the methods detectBreakpoints and mergeBreakpoints. Since only one plot is made, the function will only work for objects of class Breakpoints having length one.
geneSymbols	Boolean value whether to automatically load and plot the gene symbols from the Ensembl database. Additionally, geneSymbols can be a vector of two strings for an own annotation.
plotMut	Boolean value whether to mark deletions, insertions and mismatches.
plotBasePair	S
	Optionally, plotChimericReads displays all base pairs in a given region around the breakpoint (see maxBasePairs).
maxBasePairs	The maximum number of base pairs to be plotted. Only used in conjunction with plotBasePairs=TRUE.
legend	A logical value (TRUE/FALSE) whether to plot a legend that explains the colouration of the insertions, deletions, mismatches and breakpoints.
title	A title for the plot.
col	A vector of four colours to draw insertions, deletions, mismatches and breakpoints. In this order, the default colours are "red", "green", "black" and "orange" (use colours () to see a list of possible values).

Details

This method is intended to be run after the pipeline for structural variant detection. Therefore, see the methods filterChimericReads, detectBreakpoints and mergeBreakpoints to correctly preprocess your alignment before running plotChimericReads.

Note

It is recommended to first create and resize the output device (e.g. the plotting window or a pdf file) before plotting. For example, on Unix systems you may try X11 (width=w, height=h) or pdf (file="plotChimericReads.pdf", width=w, height=h) for some window width w (e.g. w=12) and window height h (e.g. h=6).

plot Variants 47

Author(s)

Christoph Bartenhagen

See Also

Breakpoints-class, detectBreakpoints, mergeBreakpoints

Examples

```
# load breakpoint data containing twelve chimeric reads describing an inversion in chromo
data("breakpoints")
breakpoints

# standard plot
# (only arrangement of reads plotted; breakpoints in orange, deletions
# in red, insertions in green and mismatches in black by default)
plotChimericReads(breakpoints)

# plot base pairs in the breakpoint region (+/- 32bp)
## Not run: plotChimericReads(breakpoints, plotBasePairs=TRUE, maxBasePairs=32)

# use custom colours and display a legend:
# deletions="brown", insertions="blue", mismatches="yellow", breakpoints="gray"
plotChimericReads(breakpoints, col=c("brown", "blue", "yellow", "gray"), legend=TRUE)
```

plotVariants

Plots variant positions

Description

This function illustrates the positions and types of all variants within a given transcript.

Usage

```
plotVariants(varData, transcript, legend=TRUE, regions=c(), regLabel="(region la
```

Arguments

varData An instance of class annotated Variants you get by calling annotate Variants.

If no such argument is supplied the data will be read from the Ensembl database

automatically for all variants.

transcript This argument can either be a string containing an Ensembl transcript-id or a

data frame containing the associated Ensembl transcript information. In the latter case, the data frame requires the columns "ensembl_transcript_id", "rank", "cds_start", "cds_end" and "cds_length". If only the Ensembl transcript-id is

passed, the function will return an appropriate data frame.

legend A logical value (TRUE/FALSE) whether to plot a legend that explains the type

of mutation.

regions A vector with transcript positions (start and end for every region) to be high-

lighted in the plot. It must have the form regions=c(start1, end1,

start2, end2, start3, end3 ...).

48 plot Variants

regLabel	If legend=TRUE and regions were specified, it is recommended to give a short label that explains the highlighted region(s).
col	A vector specifying four colours for missense point mutations, silent point mutations, nonsense point mutations and deletions (in this order). By default, these mutations are drawn as grey, white, black arrows and grey diamonds respectively.
title	A title for the plot.

Details

The plot shows only coding regions and distinguishes between missense, silent and nonsense point mutations and deletions.

Value

If an Ensembl transcript-id is passed as transcript argument, the function will return a data frame containing the Ensembl transcript information required for plotting ("ensembl_transcript_id", "rank", "cds_start", "cds_end" and "cds_length"). This data frame may prove useful for future plots with the same transcript.

If such a data frame is passed directly, the function returns nothing.

Author(s)

Christoph Bartenhagen

See Also

```
annotateVariants.
```

```
# one missense, one nonsense point mutation and one deletion
variants = data.frame(
    row.names=c("missense", "deletion", "nonsense"),
    start=c(106157528, 106157635, 106193892),
    end=c(106157528, 106157635, 106193892),
    chromosome=c("4", "4", "4"),
    strand=c("+", "+", "+"),
    seqRef=c("A", "G", "C"),
    seqMut=c("G", "-", "T"),
    seqSur=c("TACAGAA", "TAAGCAG", "CGGCGAA"),
    stringsAsFactors=FALSE)

# annotate variants with affected genes, exons and codons (may take a minute to finish)
## Not run: varAnnot = annotateVariants(variants)

# plot variants vor transcript "ENST00000380013" and mark region within [700, 1000]
## Not run: plotVariants(varAnnot, transcript="ENST00000380013", legend=TRUE,regions=c(70)
```

```
plotVariationFrequency
```

Create an AVA style variation frequency plot

Description

This method creates a plot similar to the variation frequency plot in Roche's GS Amplicon Variant Analyzer. The plot shows the reference sequence along the x-axis and indicates variants as bars at the appropriate positions. The height of the bars corresponds to the percentage of reads carrying the variant. A second y-axis indicates the absolute number of reads covering the variant.

Usage

```
plotVariationFrequency(object, plotRange, ...)
```

Arguments

object	A character pointing to an Amplicon Variant Analyser Global Alignmnet export file.
plotRange	A two dimensional numeric vector giving the start and end base of the reference sequence that should be plotted.
•••	Arguments passed to other plotting methods. Especially, argument col: Optional character vector of length 7 specifying the bars' colors indicating different base substitutions or deletions. See details. And argument sequenceCex: Optional numeric value specifying the size of the reference sequence's bases.

Details

The text file used as imput must have the format generated by the AVA export function. Such a file can be generated using the export button in the Global Alignment view of the AVA software. The col argument specifies the colours used for different bases and deletions. The following listing gives the meaning of the i-th position of the col vector (default values in braces):

- 1. A (green)
- 2. C (blue)
- 3. G (black)
- 4. T (red)
- 5. N (purple)
- 6. deletion (gray)

Author(s)

Hans-Ulrich Klein

```
## Not run:
file = system.file("extdata", "AVAVarFreqExport", "AVAVarFreqExport.xls", package="R453Pl
plotVariationFrequency(file, plotRange=c(50, 150))
## End(Not run)
```

50 readLengthHist

```
positionQualityBoxplot
```

Boxplot Of The Quality For Each Position

Description

Creates a boxplot of the quality scores over all sequences at each position.

Usage

```
positionQualityBoxplot(object, range, binsize=10,
    xlab=paste("Read position in bp (Bin size: ", binsize, "bp)", sep=""), ylab=
    col="firebrick1", ...)
```

Arguments

object	An object of class QualityScaledDNAStringSet, ShortReadQ or SFFContainer.
range	A numeric vector of length one or two. If length one only bases from the first until this position are plotted. If two all bases between these two positions are plotted.
binsize	Number of positions to summarize in one box in the plot.
xlab	The X axis label.
ylab	The Y axis label.
col	The plotting color.
	Arguments to be passed to methods, such as graphical parameters (see 'par').

Author(s)

Christian Ruckert

```
readLengthHist Histogram Of The Read Lengths
```

Description

This function plots a histogram of the read lengths.

Usage

```
readLengthHist(object, cutoff=0.99, xlab="Read length", ylab="Number of sequential", breaks=100, ...)
```

readLengthStats 51

Arguments

object	An object of class DNAStringSet, ShortRead or SFFContainer.
cutoff	Reads longer than the cutoff-percent quantile are omitted from the plot.
xlab	The X axis label.
ylab	The Y axis label.
col	The plotting color.
breaks	The number of breaks in the histogram (see 'hist').
	Arguments to be passed to methods, such as graphical parameters (see 'par').

Author(s)

Christian Ruckert

readLengthStats Statistics For The Read Lengths

Description

This function returns the mean, median, minimum, maximum and standard deviation of the read lengths over a set of sequences.

Usage

readLengthStats(object)

Arguments

object An object of class DNAStringSet, ShortRead or SFFContainer.

Value

A vector with five entries: mean, median, min, max and sd.

Author(s)

Christian Ruckert

52 readsOnTarget

readSff

Function To Read In Roche's .sff Files

Description

This function reads in files in Roche's Standard Flowgram Format (SFF) and store the contents in an SFFContainer-class object.

Usage

```
readSFF (files)
```

Arguments

files

The name of the .sff file to read in or a character vector of multiple file names or the name of a directory containing .sff files.

Value

An object or a list of objects of class SFFContainer storing all the information from the .sff file(s).

Author(s)

Christian Ruckert

See Also

SFFContainer.

Examples

```
## Not run: sffContainer = readSFF("test.sff")
```

readsOnTarget

Check for each read whether it aligns within the given region.

Description

This methods checks (approximately) whether the given reads align within a given region.

Usage

```
readsOnTarget(alnReads, targetRegion)
```

Arguments

alnReads A list as returned by scanBam storing aligned reads.

targetRegion The target region as a RangesList. The chromosome names must fit to the chromosome names used in the alignment information of the given reads.

referenceSequences 53

Details

The detailed alignment information given by the CIGAR strings in .bam files are ignored by the function. Instead, it is assumed that the whole read alignes to the reference without indels. This is often not true for longer read (e.g. generated with Roche 454 Sequencing), but saves computation time. Hence, this method is useful to approximate the number of reads that align in the target region of a targeted sequencing experiment.

Value

A list with one logical vector for each list entry in alnReads. The logical vector indicates for each read whether it overlaps with at least one base from any target region or not.

Author(s)

Hans-Ulrich Klein

See Also

scanBam

Examples

```
library(Rsamtools)
bamFile = system.file("extdata", "StructuralVariantDetection", "bam", "N01.bam", package=
bam = scanBam(bamFile)
region = RangesList("11"=IRanges(start=118307205, end=118395936))
targetReads = readsOnTarget(bam, region)
sum(targetReads[[1]])
```

referenceSequences Access the reference sequences of an AVASet

Description

This function give access to a slot of an instance of the AVASet storing information about all reference sequences of the amplicons.

Usage

```
referenceSequences (object)
```

Arguments

```
object An link (AVASet-class) object.
```

Value

The data is stored in an object of class AlignedRead and thus gives information about all reference sequences and their position on a chromosome (if alignShortReads has been called before).

54 removeLinker

Author(s)

Christoph Bartenhagen

See Also

```
alignShortReads
```

Examples

```
# load an AVA dataset containing 6 samples, 4 amplicons and 259 variants
data(avaSetExample)
```

referenceSequences (avaSetExample)

removeLinker

Remove linker sequences located at the start of short reads

Description

If linkers are attached during sample preparation, it may be useful to remove the linkers' sequences after sequencing. This method finds and removes linker sequences that are located at the start of the given reads.

Usage

```
## S4 method for signature 'XStringSet, DNAString, logical, numeric, numeric'
removeLinker(reads, linker, removeReadsWithoutLinker, minOverlap, penalty)
```

Arguments

reads A DNAStringSet instance that contains reads possibly having linkers at their

start site

linker A DNAString instance with the linker's sequence

removeReadsWithoutLinker

Whether reads without linkers should be removed. Default is FALSE

minOverlap The minimal score that must be achived when aligning the linker. Default is

length(linker)/2

penalty The penalty for substitutions or indels. Default is -2

Details

The best alignment of the linker within the start (length of linker + 5) of each given sequence is computed. The followong scoring schema is used: Each matching bases scores +1. Each substitution or indel scores the given penalty argument (default: penalty=-2). There are no penalties for gaps and the end of the linker (overlap). An alignment is considered as match, if the scores is larger of equal to minOverlap (default: minOverlap=round(length(linker)/2)). In cases of a successful match, the subsequence from position 1 until the end of the linker's alignment is removed.

Value

removeLinker returns a DNAStringSet with trimmed reads.

Author(s)

Hans-Ulrich Klein

See Also

```
sequenceCaptureLinkers, DNAStringSet, pairwiseAlignment
```

Examples

sequenceCaptureLinkers

Retrieve NimbleGen's sequence capture linkers

Description

This method returns the NimbleGen's linker sequences used with their sequence capture arrays. See pp.29-30 in the NimbleGen Arrays User's Guide.

Usage

Arguments

name

Character vector with linker sequences' names

Details

If the argument name is omitted, both linker sequences are returned.

Value

 ${\tt sequenceCaptureLinkers} \ a \ {\tt DNAStringSet} \ with \ the \ requested \ linker \ sequences.$

Author(s)

Hans-Ulrich Klein

56 setVariantFilter

See Also

```
removeLinker
```

Examples

```
sequenceCaptureLinkers()
```

```
sequenceQualityHist
```

A Histogram Of The Sequence Qualities

Description

This function creates a histogram of the mean qualities of the sequences.

Usage

```
sequenceQualityHist(object, xlab="Mean of quality scores per sequence",
  ylab="Number of sequences", col="firebrick1", ...)
```

Arguments

object	An object of class QualityScaledDNAStringSet, ShortReadQ or SFFContainer.
xlab	The X axis label.
ylab	The Y axis label.
col	The plotting color.
	Arguments to be passed to methods, such as graphical parameters (see 'par').

Author(s)

Christian Ruckert

```
setVariantFilter Filters output of variant information
```

Description

This functions sets the filter to display only those variants, whose amplicon coverage (in percent) in forward and reverse direction in at least one sample is higher than a given value. The coverage is defined as the percentual amount of reads that cover a variant.

Usage

```
setVariantFilter(object, filter=0)
```

setVariantFilter 57

Arguments

object	An instance of an link {AVASet-class} or MapperSet-class.
filter	A filter value between 0 and 1. If two values are given in a vector, the variants are filtered according to the forward (first value) and reverse direction (second value) separately. In this case, a variant has to meet both requirements.

Details

Setting the filter affects the assayData and the featureData of the variant slot. See also getVariantPercentages for further details.

Value

```
setVariantFilter returns the given link {AVASet-class}/link {MapperSet-class} instance with an updated filter value.
```

Author(s)

Christoph Bartenhagen

See Also

```
link{AVASet-class}, link{MapperSet-class}, getVariantPercentages.
```

```
# load an AVA dataset containing 6 samples, 4 amplicons and 259 variants
data(avaSetExample)
avaSetExample

# use only those variants that are covered by at least 10% of all reads in one sample in
avaSetExample = setVariantFilter(avaSetExample, filter=0.1)
avaSetExample

# use only those variants that are covered by at least 0.1% of all reads in one sample in
# and by at least 0% in reverse direction (259 -> 6 variants)
avaSetExample = setVariantFilter(avaSetExample, filter=c(0.1, 0))
avaSetExample

# reset filter values to zero
avaSetExample = setVariantFilter(avaSetExample, filter=0)
# or simply
avaSetExample = setVariantFilter(avaSetExample)
```

58 transcriptdf

Write A SFFContainer Object To A FASTQ File

Description

This function takes a SFFContainer object and writes it to a file in FASTQ format.

Usage

```
sff2fastq(x, outdir, fname)
```

Arguments

x An object of class SFFContainer.

outdir The directory where the file should be stored, defaults to the current working

directory.

fname The name of the file to write. Defaults to the filename slot of the SFFContainer,

with .sff substituted with .fastq.

Author(s)

Christian Ruckert

transcriptdf

Ensembl transcript information for variant plotting

Description

This example contains Ensembl transcript information (here for transcript 'ENST00000380013') as created, used and returned by the function link{plotVariants}. It mainly holds locations of coding areas within the transcript that will then be highlighted in the variant plot.

Usage

```
data(transcriptdf)
```

Format

A data frame with 9 observations on the following 5 variables.

```
ensembl_transcript_id Ensembl id for the transcript
rank exon rank
cds_start start of a codon
cds_end end of a codon
cds_length length of a codon
```

```
data(transcriptdf)
```

Index

*Topic TiTv,Transition,Transversion	homopolymerHist, 40
calculateTiTv, 22	nucleotideCharts,44
*Topic Variation Frequency Coverage	positionQualityBoxplot,50
plotVariationFrequency, 49	readLengthHist, 50
*Topic alignShortReads	readLengthStats, 51
alignShortReads, 15	sequenceQualityHist,56
*Topic amplicon coverage	sff2fastq, 58
plotAmpliconCoverage, 45	*Topic plotChimericReads,
*Topic annotateVariants	detectBreakpoints,
annotateVariants, 16	mergeBreakpoints,
*Topic classes	Breakpoints
AnnotatedVariants-class, 5	plotChimericReads, 46
AVASet-class, 1	*Topic readsOnTarget
Breakpoints-class, 6	readsOnTarget, 52
MapperSet-class,8	*Topic removeLinker
SFFContainer-class, 11	removeLinker, 54
SFFRead-class, 13	[, AVASet, ANY, ANY-method
*Topic datasets	(AVASet-class), 1
avaSetExample, 18	[,Breakpoints,ANY,ANY-method
avaSetFiltered, 19	(Breakpoints-class), 6
avaSetFiltered_annot, 19	[,SFFContainer,ANY,ANY-method
breakpoints, 22	(SFFContainer-class), 11
captureArray, 23	112 1/2222 / / 1 11
mapperSetExample, 42	addRead(SFFContainer-class), 11
transcriptdf, 58	addRead, SFFContainer, SFFRead-method
*Topic demultiplex	(SFFContainer-class), 11
demultiplexReads, 27	AlignedRead, 16, 30
genomeSequencerMIDs, 37	AlignedRead-class, 6
*Topic filterChimericReads	alignedReadsC1 (Breakpoints-class), 6
filterChimericReads, 33	alignedReadsC1,Breakpoints-method
*Topic linker	(Breakpoints-class), 6
sequenceCaptureLinkers,55	alignedReadsC1<-
*Topic methods	(Breakpoints-class), 6
baseFrequency, 20	alignedReadsC1<-,Breakpoints,list-method
baseQualityHist, 21	(Breakpoints-class), 6
baseQualityStats, 21	alignedReadsC2
complexity.dust, 24	(Breakpoints-class), 6
complexity.entropy, 25	alignedReadsC2, Breakpoints-method
dinucleotideOddsRatio, 31	(Breakpoints-class), 6
flowgramBarplot, 35	alignedReadsC2<-
gcContent, 35	(Breakpoints-class), 6
gcContentHist, 36	alignedReadsC2<-,Breakpoints,list-method
gcPerPosition, 36	(Breakpoints-class), 6

```
alignShortReads, 2, 4, 15, 53, 54
                                          AVASet, 3, 9, 16, 45
alignShortReads, AVASet, BSgenome, chara AVASetoginaalaquethouethod (AVASet),
       (alignShortReads), 15
alignShortReads, AVASet, BSgenome, chara ctiAscatiascs, nleth, d8, 32, 33, 38, 39,
                                                  41
       (alignShortReads), 15
alignShortReads, AVASet, BSgenome, missinAy ASegical and thod
       (alignShortReads), 15
                                          avaSetExample, 18
alignShortReads, AVASet, BSgenome, missing Anstring anech 10
                                          avaSetFiltered_annot, 19
       (alignShortReads), 15
alignShortReads, AVASet, DNAStringSet, character-method
                                          baseFrequency, 20
       (AVASet-class), 1
alignShortReads,DNAStringSet,BSgenome,baseFrequency,DNAStringSet-method
                                                  (baseFrequency), 20
       (alignShortReads), 15
alignShortReads, DNAStringSet, BSgenome, baseFrequency, SFFContainer-method
                                                  (baseFrequency), 20
       (alignShortReads), 15
alignShortReads, DNAStringSet, BSgenome, missing, logical-method
                                                  (baseFrequency), 20
       (alignShortReads), 15
alignShortReads, DNAStringSet, BSgenome, missing, mrssing method baseQualityHist, QualityScaledDNAStringSet-method
       (alignShortReads), 15
                                                  (baseQualityHist), 21
alphabetFrequency, 20
                                          baseQualityHist,SFFContainer-method
AnnotatedDataFrame, 33
                                                  (baseQualityHist), 21
annotatedVariants
                                          baseQualityHist,ShortReadQ-method
       (AnnotatedVariants-class).
                                                  (baseQualityHist), 21
\label{eq:annotatedVariants} \begin{array}{c} \text{baseQualityStats, } \mathbf{21} \\ \text{annotatedVariants-method} \end{array}
                                           baseQualityStats,QualityScaledDNAStringSet-met
       (AnnotatedVariants-class),
                                                  (baseQualityStats), 21
                                          baseQualityStats,SFFContainer-method
AnnotatedVariants-class, 17
                                                  (baseQualityStats), 21
AnnotatedVariants-class, 5
                                          baseQualityStats,ShortReadQ-method
annotatedVariants<-, AnnotatedVariants, list-method (baseQualityStats), 21
       (AnnotatedVariants-class),
                                          breakpoints, 22
                                          Breakpoints-class, 34, 44, 47
annotateVariants, 2, 5, 9, 16, 41, 42, 47,
                                          Breakpoints-class, 6
annotateVariants, AVASet, missing-methodalculateTiTv, 22
       (annotateVariants), 16
                                          calculateTiTv, AVASet-method
annotateVariants, AVASet-method
                                                  (calculateTiTv), 22
       (AVASet-class), 1
                                          calculateTiTv, MapperSet-method
annotateVariants, data.frame, missing-method (calculateTiTv), 22
       (annotateVariants), 16
                                          captureArray, 23
annotateVariants, MapperSet, BSgenome-methodAdapterLeft
       (annotateVariants), 16
                                                  (SFFContainer-class), 11
annotateVariants, MapperSet, missing-methodAdapterLeft, SFFContainer-method
       (annotateVariants), 16
                                                  (SFFContainer-class), 11
annotateVariants, MapperSet-method
                                          clipAdapterLeft, SFFRead-method
       (MapperSet-class), 8
                                                  (SFFRead-class), 13
assayDataAmp, 17, 32, 33
                                          clipAdapterLeft<-
assayDataAmp, AVASet-method
                                                  (SFFContainer-class), 11
       (AVASet-class), 1
                                          clipAdapterLeft<-, SFFContainer, numeric-method</pre>
assayDataAmp<- (AVASet-class), 1
                                                  (SFFContainer-class), 11
assayDataAmp<-, AVASet, AssayData-methoctlipAdapterLeft<-, SFFRead, numeric-method
       (AVASet-class), 1
                                                  (SFFRead-class), 13
```

commonBpsC1 (Breakpoints-class), 6

clipAdapterRight

(SFFContainer-class), 11	commonBpsC1,Breakpoints-method
clipAdapterRight,SFFContainer-method	(Breakpoints-class), 6
(SFFContainer-class), 11	commonBpsC1<-
clipAdapterRight,SFFRead-method	(Breakpoints-class), 6
(SFFRead-class), 13	commonBpsC1<-,Breakpoints,list-method
clipAdapterRight<-	(Breakpoints-class), 6
(SFFContainer-class), 11	commonBpsC2 ($Breakpoints-class$), 6
clipAdapterRight<-,SFFContainer,numer	pommonRosC2,Breakpoints-method
(SFFContainer-class), 11	(Breakpoints-class), 6
clipAdapterRight<-,SFFRead,numeric-me	tangunan Bps C2 <-
(SFFRead-class), 13	(Breakpoints-class), 6
clipQualityLeft	commonBpsC2<-,Breakpoints,list-method
(SFFContainer-class), 11	(Breakpoints-class), 6
clipQualityLeft,SFFContainer-method	complexity.dust, 24
(SFFContainer-class), 11	complexity.dust,DNAStringSet-method
clipQualityLeft,SFFRead-method	(complexity.dust), 24
(SFFRead-class), 13	complexity.dust,SFFContainer-method
clipQualityLeft<-	(complexity.dust), 24
(SFFContainer-class).11	complexity.dust,ShortRead-method
clipQualityLeft<-,SFFContainer,numeri	c-metho(complexity.dust), 24
(CEEContainon alaga) 11	COMPLEXITY, ENTROPY, 23
clipOualityLeft <sffread.numeric-met< td=""><td>complexity.entropy, DNAStringSet-method</td></sffread.numeric-met<>	complexity.entropy, DNAStringSet-method
(SFFRead-class), 13	(complexity.entropy), 25
clipQualityRight	complexity.entropy,SFFContainer-method
(SFFContainer-class), 11	(complexity.entropy), 25
clipQualityRight, SFFContainer-method	complexity.entropy,ShortRead-method
(SFFContainer-class), 11	(complexity.entropy), 25
clipQualityRight,SFFRead-method	convertCigar, 26
(SFFRead-class), 13	coverageOnTarget, 26
clipQualityRight<-	<pre>coverageOnTarget,list,RangesList-method</pre>
(SFFContainer-class), 11	(coverageOnTarget), 26
	cia-mothod
<pre>clipQualityRight<-,SFFContainer,numer</pre>	
	DataFrame, 29
<pre>clipQualityRight<-,SFFRead,numeric-me (SFFRead-class),13</pre>	
commonAlignC1	demultiplexReads, XStringSet, XStringSet, missing
_	(demultiplexReads), 27
(Breakpoints-class), 6 commonAlignC1,Breakpoints-method	demultiplexReads, XStringSet, XStringSet, missing
(Breakpoints-class), 6	(demultiplexReads), 27
	demultiplexReads, XStringSet, XStringSet, numeric
commonAlignC1<-	(demultiplexReads), 27
(Breakpoints-class), 6	demultiplexReads, XStringSet, XStringSet, numeric
commonAlignC1<-,Breakpoints,list-meth	(1
(Breakpoints-class), 6	detectBreakpoints, 6, 8, 28, 34, 43, 44,
commonAlignC2	46, 47
(Breakpoints-class), 6	detectBreakpoints, list-method
commonAlignC2, Breakpoints-method	(detectBreakpoints), 28
(Breakpoints-class), 6	dinucleotideOddsRatio,31
commonAlignC2<-	dinucleotideOddsRatio,DNAStringSet-method
(Breakpoints-class), 6	(dinucleotideOddsRatio), 31
	nochinucleotideOddsRatio,SFFContainer-method
(Breakpoints-class),6	(dinucleotideOddsRatio), 31

dinucleotideOddsRatio,ShortRead-metho	oflowgram<-,SFFRead,numeric-method
(dinucleotideOddsRatio),31	(SFFRead-class), 13
DNAStringSet, 16, 20, 24, 25, 28, 31,	flowgramBarplot, 35
35–38, 44, 51, 55	flowgramBarplot, SFFRead-method
	(flowgramBarplot), 35
eSet, 2, 9	flowgramFormat
extendedCIGARToList, 26	(SFFContainer-class), 11
extendedCIGARToList	flowgramFormat, SFFContainer-method
(convertCigar), 26	(SFFContainer-class), 11
	flowgramFormat, SFFRead-method
fData, 41	(SFFRead-class), 13
fDataAmp, 18, 32, 33, 38	flowgramFormat<-
fDataAmp, AVASet-method	(SFFContainer-class), 11
(AVASet-class), 1	<pre>flowgramFormat<-,SFFContainer,numeric-method</pre>
featureDataAmp, 18, 32, 32	(SFFContainer-class), 11
featureDataAmp, AVASet-method	flowgramFormat<-,SFFRead,numeric-method
(AVASet-class), 1	(SFFRead-class), 13
featureDataAmp<- (AVASet-class) 1	flowgrams(SFFContainer-class),11
featureDataAmp<-, AVASet, AnnotatedData	Flamermethod
(AVASet-class), 1	(SFFContainer-class), 11
filterChimericReads, 8, 28-30, 33, 46	flowgrams<-(SFFContainer-class),
<pre>filterChimericReads, list, missing, DNAS</pre>	string,mlssing,missing-method flowgrams<-,SFFContainer,list-method
filterChimericReads, list, missing, DNAS (filterChimericReads), 33	tring, numeric, numeric hats) old flowIndexes(SFFContainer-class),
filterChimericReads, list, missing, miss	ing, missing, missing-method
<pre>(filterChimericReads), 33 filterChimericReads, list, missing, miss</pre>	
(filterChimericReads),33	llowindexes, Sff Read-Method
<pre>filterChimericReads, list, RangesList, D</pre>	NAStrin GSATESATESATES ing-method flowIndexes<-
filterChimericReads, list, RangesList, D	NAString, Franceria, namerias method flowIndexes<-, SFFContainer, list-method
<pre>(filterChimericReads), 33 filterChimericReads, list, RangesList, m</pre>	nissing, MFFFG99t, minesinglamethold
(filterChimericReads), 33	flowIndexes<-,SFFRead,numeric-method
<pre>filterChimericReads, list, RangesList, m</pre>	
flowChars(SFFContainer-class),11	gcContent, 35
flowChars,SFFContainer-method	gcContent,DNAStringSet-method
(SFFContainer-class), 11	(gcContent), 35
flowChars,SFFRead-method	gcContent,SFFContainer-method
(SFFRead-class), 13	(gcContent), 35
flowChars<-(SFFContainer-class),	gcContent,ShortRead-method
11	(gcContent), 35
<pre>flowChars<-,SFFContainer,character-me</pre>	topoContentHist,36
(SFFContainer-class), 11	gcContentHist,DNAStringSet-method
<pre>flowChars<-, SFFRead, character-method</pre>	(gcContentHist), 36
(SFFRead-class), 13	gcContentHist,SFFContainer-method
flowgram (SFFRead-class), 13	(gcContentHist), 36
flowgram, SFFRead-method	gcContentHist,ShortRead-method
(SFFRead-class), 13	(gcContentHist), 36
flowgram<-(SFFRead-class), 13	gcPerPosition, 36

gcPerPosition,DNAStringSet-method	listToExtendedCIGAR
(gcPerPosition), 36	(convertCigar), 26
gcPerPosition,SFFContainer-method	
(gcPerPosition), 36	MapperSet, 9, 16
gcPerPosition,ShortRead-method	MapperSet, character-method
(gcPerPosition), 36	(MapperSet), 9
genomeSequencerMIDs, 28, 37	MapperSet-class, 2, 4, 17, 39, 41, 57
genomeSequencerMIDs, character-method	
(genomeSequencerMIDs), 37	mapperSetExample, 42
genomeSequencerMIDs, missing-method	matchPDict, 16
(genomeSequencerMIDs), 37	mergeBreakpoints, 6, 8, 29, 30, 34, 43,
getAlignedReads, 38	46, 47
getAlignedReads, AVASet-method	mergeBreakpoints, Breakpoints, missing, list-met
(getAlignedReads), 38	(Breakpoints-class), 6
getAminoAbbr, 39	mergeBreakpoints, Breakpoints, missing, missing-
getRead, 13	(Breakpoints-class), 6
getRead(<i>SFFContainer-class</i>),11	mergeBreakpoints, Breakpoints, numeric, missing-
getRead, SFFContainer, character-method	(Bleakpoints-class), 0
(SFFContainer-class), 11	name (SFFContainer-class), 11
getReadStatus (MapperSet-class), 8	name, SFFContainer-method
getReadStatus, MapperSet-method	(SFFContainer-class), 11
(MapperSet-class), 8	
getVariantPercentages, 2, 9, 39, 57	name, SFFRead-method
getVariantPercentages, AVASet-method	(SFFRead-class), 13
(AVASet-class), 1	name<-(SFFContainer-class), 11
	omame<-, SFFContainer, character-method
(MapperSet-class),8	(SFFContainer-class), 11
homonolymorujat 40	name<-, SFFRead, character-method
homopolymerHist, 40	(SFFRead-class), 13
homopolymerHist, SFFRead-method	names, Annotated Variants-method
(homopolymerHist), 40	(AnnotatedVariants-class),
htmlReport, 2, 5, 9, 17, 41	5
htmlReport, AVASet-method	names, Breakpoints-method
(AVASet-class), 1	(Breakpoints-class), 6
htmlReport, MapperSet-method	names<-, Annotated Variants, character-method
(MapperSet-class),8	(AnnotatedVariants-class),
IRanges, 29	5
11.dilgC3, 27	names<-,Breakpoints,ANY-method
keySequence(SFFContainer-class),	(Breakpoints-class), 6
11	nucleotideCharts,44
keySequence, SFFContainer-method	nucleotideCharts,DNAStringSet-method
(SFFContainer-class), 11	(nucleotideCharts),44
keySequence, SFFRead-method	nucleotideCharts,SFFContainer-method
(SFFRead-class), 13	(nucleotideCharts),44
keySequence<-	nucleotideCharts, ShortRead-method
(SFFContainer-class), 11	(nucleotideCharts),44
keySequence<-, SFFContainer, character-	-method
(SEEContainer-class) 11	PairwiseAlignedFixedSubject, 30
keySequence<-,SFFRead,character-metho	_PairwiseAlignedFixedSubject-
(SFFRead-class), 13	class,
(SEFFICAU CLASS), 13	6
length, Breakpoints-method	pairwiseAlignment, 55
(Breakpoints-class), 6	plotAmpliconCoverage, 45
	=

```
plotAmpliconCoverage, AVASet, character, respectives that the Strates, DNAStringSet-method
       (plotAmpliconCoverage), 45
                                                 (readLengthStats), 51
plotAmpliconCoverage, AVASet, character, meiasticing time that still states, SFFContainer-method
       (plotAmpliconCoverage), 45
                                                 (readLengthStats), 51
plotAmpliconCoverage, AVASet, missing, logical engeth States, ShortRead-method
       (plotAmpliconCoverage), 45
                                                 (readLengthStats), 51
plotAmpliconCoverage, AVASet, missing, missing SNFCood ainer-class), 11
       (plotAmpliconCoverage), 45
                                         reads, SFFContainer-method
plotChimericReads, 8, 30, 44, 46
                                                 (SFFContainer-class), 11
plotChimericReads, Breakpoints-method reads<-(SFFContainer-class), 11
       (Breakpoints-class), 6
                                         reads<-, SFFContainer, QualityScaledDNAStringSet
plotVariants, 47
                                                 (SFFContainer-class), 11
plotVariants, AnnotatedVariants, charactreads th, 44, 13, 15
       (plotVariants), 47
                                         readSFF (readSff), 52
plotVariants, Annotated Variants, data.framegynethod
       (plotVariants), 47
                                         readsOnTarget, 52
plotVariationFrequency, 49
                                         readsOnTarget,list,RangesList-method
plotVariationFrequency, character, numeric-methedadsOnTarget), 52
       (plotVariationFrequency),
                                         referenceSequences, 53
                                         referenceSequences, AVASet-method
positionQualityBoxplot, 50
                                                 (AVASet-class), 1
positionQualityBoxplot,QualityScaledDNAStringSet-quethed<-
       (positionQualityBoxplot),
                                                 (AVASet-class), 1
                                         referenceSequences<-, AVASet, AlignedRead-method
positionQualityBoxplot,SFFContainer-method
                                                (AVASet-class), 1
       (positionQualityBoxplot),
                                         removeLinker, 54, 56
                                         removeLinker, XStringSet, DNAString, logical, nume
positionQualityBoxplot,ShortReadQ-method
                                                 (removeLinker), 54
       (positionQualityBoxplot),
                                          removeLinker, XStringSet, DNAString, missing, miss
       50
                                                 (removeLinker), 54
quality, SFFRead-method
                                         scanBam, 27, 28, 33, 34, 53
       (SFFRead-class), 13
                                         seqsC1 (Breakpoints-class), 6
quality<- (SFFRead-class), 13
                                         seqsC1, Breakpoints-method
quality<-, SFFRead, BString-method
                                                 (Breakpoints-class), 6
       (SFFRead-class), 13
                                         seqsC1<-(Breakpoints-class), 6
QualityScaledDNAStringSet, 21, 50,
                                         seqsC1<-,Breakpoints,list-method</pre>
       56
                                                 (Breakpoints-class), 6
                                         seqsC2 (Breakpoints-class), 6
read (SFFRead-class), 13
                                         seqsC2, Breakpoints-method
read, SFFRead-method
                                                 (Breakpoints-class), 6
       (SFFRead-class), 13
                                         seqsC2<-(Breakpoints-class), 6</pre>
read<-(SFFRead-class), 13
read<-, SFFRead, QualityScaledDNAStringSetgan2tmo&reakpoints, list-method
                                                 (Breakpoints-class), 6
       (SFFRead-class), 13
                                         sequenceCaptureLinkers, 34, 55, 55
readLengthHist, 50
                                          sequenceCaptureLinkers, character-method
readLengthHist,DNAStringSet-method
                                                 (sequenceCaptureLinkers),
       (readLengthHist), 50
                                                 55
readLengthHist, SFFContainer-method
       (readLengthHist), 50
                                          sequenceCaptureLinkers, missing-method
readLengthHist,ShortRead-method
                                                 (sequenceCaptureLinkers),
       (readLengthHist), 50
readLengthStats, 51
                                          sequenceQualityHist, 56
```

```
sequenceQualityHist,QualityScaledDNAStringSet-method
       (sequenceQualityHist), 56
sequenceQualityHist,SFFContainer-method
       (sequenceQualityHist), 56
sequenceQualityHist,ShortReadQ-method
       (sequenceQualityHist), 56
setVariantFilter, 1, 2, 8, 9, 19, 39, 56
setVariantFilter, AVASet-method
       (AVASet-class), 1
setVariantFilter,MapperSet-method
       (MapperSet-class), 8
sff2fastq, 58
sff2fastq, SFFContainer-method
       (sff2fastq), 58
SFFContainer, 12, 13, 15, 20, 21, 24, 25,
       31, 35–37, 44, 50–52, 56, 58
SFFContainer
       (SFFContainer-class), 11
SFFContainer-class, 52
SFFContainer-class, 11
SFFRead, 12, 13, 35, 40
SFFRead (SFFRead-class), 13
SFFRead-class, 13
ShortRead, 20, 24, 25, 31, 35-37, 44, 51
ShortReadQ, 21, 50, 56
transcriptdf, 58
vector, 51
Versioned, 2, 9
VersionedBiobase, 2, 9
```