Package 'PeacoQC'

October 15, 2023

Title Peak-based selection of high quality cytometry data

Version 1.10.0

Description

This is a package that includes pre-processing and quality control functions that can remove margin events, compensate and transform the data and that will use PeacoQCSignalStability for quality control. This last function will first detect peaks in each channel of the flowframe. It will remove anomalies based on the IsolationTree function and the MAD outlier detection method. This package can be used for both flow- and mass cytometry data.

```
Encoding UTF-8
License GPL (>=3)
LazyData true
URL http://github.com/saeyslab/PeacoQC
BugReports http://github.com/saeyslab/PeacoQC/issues
Depends R (>= 4.0)
Imports circlize, ComplexHeatmap, flowCore, flowWorkspace, ggplot2,
     grDevices, grid, gridExtra, methods, plyr, stats, utils
RoxygenNote 7.2.0
Suggests knitr, rmarkdown, BiocStyle
VignetteBuilder knitr
biocViews FlowCytometry, QualityControl, Preprocessing, PeakDetection
git_url https://git.bioconductor.org/packages/PeacoQC
git_branch RELEASE_3_17
git_last_commit 4a9e75a
git_last_commit_date 2023-04-25
Date/Publication 2023-10-15
Author Annelies Emmaneel [aut, cre]
Maintainer Annelies Emmaneel <annelies.emmaneel@hotmail.com>
```

2 PeacoQC

R topics documented:

PeacoQC	 . 2
PeacoQCHeatmap	 . 4
PlotPeacoQC	 . 5
RemoveDoublets	 . 7
RemoveMargins	 . 8
	10

PeacoQC

Index

Peak-based detection of high quality cytometry data

Description

PeacoQC will determine peaks on the channels in the flowframe. Then it will remove anomalies caused by e.g. clogs, changes in speed etc. by using an IsolationTree and/or the MAD method.

Usage

```
PeacoQC(ff, channels, determine_good_cells="all",
    plot=20, save_fcs=TRUE, output_directory=".",
    name_directory="PeacoQC_results", report=TRUE,
    events_per_bin=FindEventsPerBin(remove_zeros, ff, channels,
    min_cells, max_bins, step), min_cells=150, max_bins=500, step=500,
    MAD=6, IT_limit=0.6, consecutive_bins=5, remove_zeros=FALSE,
    suffix_fcs="_QC", force_IT=150, peak_removal = (1/3),
    min_nr_bins_peakdetection = 10, time_channel_parameter = "Time",
    ...)
```

Arguments

ff

A flowframe or the location of an fcs file. Make sure that the flowframe is compensated and transformed. If it is mass cytometry data, only a transformation is necessary.

channels

Indices or names of the channels in the flowframe on which peaks have to be determined.

determine_good_cells

If set to FALSE, the algorithm will only determine peaks. If it is set to "all", the bad measurements will be filtered out based on the MAD and IT analysis. It can also be put to "MAD" or "IT" to only use one method of filtering.

plot

When PeacoQC removes more than the specified percentage, an overview plot will be made of all the selected channels and the deleted measurements. If set to TRUE, the PlotPeacoQC function is run to make an overview plot of the deleted measurements, even when nothing is removed. Default is set to 20. If an increasing or decreasing trend is found, a figure will also be made except if plot is set to FALSE.

PeacoQC 3

save_fcs If set to TRUE, the cleaned fcs file will be saved in the output_directory as: filename_QC.fcs. The _QC name can be altered with the suffix_fcs parameter. An extra column named "Original_ID" is added to this fcs file where the cells are given their original cell id. Default is TRUE.

output_directory

Directory where a new folder will be created that consists of the generated fcs files, plots and report. If set to NULL, nothing will be stored. The default folder is the working directory.

name_directory Name of folder that will be generated in output_directory. The default is "PeacoOC results".

report Overview text report that is generated after PeacoQC is run. If set to FALSE, no report will be generated. The default is TRUE.

events_per_bin Number of events that are put in one bin. Default is calculated based on the rows in ff

The minimum amount of cells (nonzero values) that should be present in one bin. Lowering this parameter can affect the robustness of the peak detection. Default is 150.

max_bins The maximum number of bins that can be used in the cleaning process. If this value is lowered, larger bins will be made. Default is 500.

The step in events_per_bin to which the parameter is reduced to. Default is 500.

MAD The MAD parameter. Default is 6. If this is increased, the algorithm becomes less strict.

The IsolationTree parameter. Default is 0.55. If this is increased, the algorithm becomes less strict.

consecutive_bins

IT_limit

If 'good' bins are located between bins that are removed, they will also be marked as 'bad'. The default is 5.

remove_zeros If this is set to TRUE, the zero values will be removed before the peak detection step. They will not be indicated as 'bad' value. This is recommended when cleaning mass cytometry data. Default is FALSE.

suffix_fcs The suffix given to the new fcs files. Default is "_QC".

force_IT If the number of determined bins is less than this number, the IT analysis will not be performed. Default is 150 bins.

peak_removal During the peak detection step, peaks are only kept if they are peak_removal percentage of the maximum height peak. Default is 1/3

 $min_nr_bins_peakdetection$

The percentage of number of bins in which the maximum number of peaks has to be present. Default is 10.

time_channel_parameter

Name of the time channel in ff if present. Default is "Time".

Options to pass on to the PlotPeacoQC function (display_cells, manual_cells, prefix)

4 PeacoQCHeatmap

Value

This function returns a list with a number of items. It will include "FinalFF" where the transformed, compensated and cleaned flowframe is stored. It also contains the starting parameters and the information necessary to give to PlotPeacoQC if the two functions are run seperatly. The Good-Cells list is also given where 'good' measurements are indicated as TRUE and the to be removed measurements as FALSE.

Examples

```
# General pipeline for preprocessing and quality control with PeacoQC
# Read in raw fcs file
fileName <- system.file("extdata", "111.fcs", package="PeacoQC")</pre>
ff <- flowCore::read.FCS(fileName)</pre>
# Define channels where the margin events should be removed
# and on which the quality control should be done
channels \leftarrow c(1, 3, 5:14, 18, 21)
ff <- RemoveMargins(ff=ff, channels=channels, output="frame")</pre>
# Compensate and transform the data
ff <- flowCore::compensate(ff, flowCore::keyword(ff)$SPILL)</pre>
ff <- flowCore::transform(ff,</pre>
                              flowCore::estimateLogicle(ff,
                              colnames(flowCore::keyword(ff)$SPILL)))
#Run PeacoOC
PeacoQC_res <- PeacoQC(ff, channels,</pre>
                         determine_good_cells="all",
                         save_fcs=TRUE)
```

PeacoQCHeatmap

Make overview heatmap of quality control analysis

Description

PeacoQCHeatmap will make a heatmap to display all the results generated by PeacoQC. It will include the percentages of measurements that are removed in total, by the IT method and by the MAD method. It will also show the parameters that were used during the quality control.

Usage

```
PeacoQCHeatmap(report_location, show_values=TRUE, show_row_names=TRUE, latest_tests=FALSE, title="PeacoQC report", ...)
```

PlotPeacoQC 5

Arguments

report_location

The path to the PeacoQC report generated by PeacoQC.

show_values

If set to TRUE, the percentages of removed values will be displayed on the heatmap. Default is TRUE.

show_row_names

If set to FALSE, the filenames will not be displayed on the heatmap. Default is TRUE.

latest_tests

If this is set to TRUE, only the latest quality control run will be displayed in the heatmap. Default is FALSE.

title

The title that should be given to the heatmap. Default is "PeacoQC_report".

Extra parameters to be given to the Heatmap function (eg. row_split)

Value

This function returns nothing but generates a heatmap that can be saved as pdf or png

Examples

PlotPeacoQC

Visualise deleted cells of PeacoQC

Description

PlotPeacoQC will generate a png file with on overview of the flow rate and the different selected channels. These will be annotated based on the measurements that were removed by PeacoQC. It is also possible to only display the quantiles and median or only the measurements without any annotation.

Usage

PlotPeacoQC

Arguments

ff	A flowframe	
channels	Indices of names of the channels in the flowframe that have to be displayed	
output_director	ry	
	Directory where the plots should be generated. Set to NULL if no plots need to be generated. The default is the working directory.	
display_cells	The number of measurements that should be displayed. (The number of dots that are displayed for every channel) The default is 2000.	
manual_cells	Give a vector (TRUE/FALSE) with annotations for each cell to compare the automated QC with. The default is NULL.	
title_FR	The title that has to be displayed above the flow rate figure. Default is NULL.	
display_peaks	If the result of PeacoQC is given, all the quality control results will be visualised. If set to TRUE: PeacoQC will be run and only the peaks will be displayed without any quality control. If set to FALSE, no peaks will be displayed and only the events will be displayed. Default is TRUE.	
prefix	The prefix that will be given to the generated png file. Default is "PeacoQC_".	
time_unit	The number of time units grouped together for visualising event rate. The default is set to 100, resulting in events per second for most flow datasets. Suggested to adapt for mass cytometry data.	
time_channel_parameter		
	Name of the time channel in ff if present. Default is "Time".	
•••	Arguments to be given to PeacoQC if display_peaks is set to TRUE.	

Value

This function returns nothing but generates a png file in the output_directory

Examples

RemoveDoublets 7

```
PlotPeacoQC(ff, channels, display_peaks=PeacoQC_res)
## Plot only the peaks (No quality control)
PlotPeacoQC(ff, channels, display_peaks=TRUE)
## Plot only the dots of the file
PlotPeacoQC(ff, channels, display_peaks=FALSE)
```

RemoveDoublets

Remove doublet events from flow cytometry data

Description

RemoveDoublets will remove doublet events from the flowframe based on two channels.

Usage

```
RemoveDoublets(ff, channel1="FSC-A", channel2="FSC-H", nmad=4, verbose=FALSE, output="frame")
```

Arguments

ff	A flowframe that contains flow cytometry data.
channel1	The first channels that will be used to determine the doublet events. Default is "FSC-A" $$
channel2	The second channels that will be used to determine the doublet events. Default is "FSC-H"
nmad	Bandwidth above the ratio allowed (cells are kept if their ratio is smaller than the median ratio + nmad times the median absolute deviation of the ratios). Default is 4.
verbose	If set to TRUE, the median ratio and width will be printed. Default is FALSE.
output	If set to "full", a list with the filtered flowframe and the indices of the doublet event is returned. If set to "frame", only the filtered flowframe is returned. The default is "frame".

Value

This function returns either a filtered flowframe when the output parameter is set to "frame" or a list containing the filtered flowframe and a TRUE/FALSE list indicating the margin events. An extra column named "Original_ID" is added to the flowframe where the cells are given their original cell id.

RemoveMargins 8

Examples

```
# Read in data
fileName <- system.file("extdata", "111.fcs", package="PeacoQC")
ff <- flowCore::read.FCS(fileName)
# Remove doublets
ff_cleaned <- RemoveDoublets(ff)</pre>
```

RemoveMargins

Remove margin events of flow cytometry data

Description

RemoveMargins will remove margin events from the flowframe based on the internal description of the fcs file.

Usage

```
RemoveMargins(ff, channels,
channel_specifications=NULL, output="frame")
```

Arguments

ff A flowframe that contains flow cytometry data.

channels The channel indices or channel names that have to be checked for margin events channel_specifications

A list of vectors with parameter specifications for certain channels. This parameter should only be used if the values in the internal parameters description is too strict or wrong for a number or all channels. This should be one list per channel with first a minRange and then a maxRange value in a vector. This list should have the channel name found back in colnames(flowCore::exprs(ff)). If a channel is not listed in this parameter, its default internal values will be used. The default of this parameter is NULL.

output

If set to "full", a list with the filtered flowframe and the indices of the margin event is returned. If set to "frame", only the filtered flowframe is returned. The default is "frame".

Value

This function returns either a filtered flowframe when the output parameter is set to "frame" or a list containing the filtered flowframe and a TRUE/FALSE list indicating the margin events. An extra column named "Original_ID" is added to the flowframe where the cells are given their original cell id.

RemoveMargins 9

Examples

Index

```
PeacoQC, 2
PeacoQCHeatmap, 4
PlotPeacoQC, 5
RemoveDoublets, 7
RemoveMargins, 8
```